



USER MANUAL

BK105 Chill Tray




*****WARNING*****

Watch for water build up underneath unit due to condenser defrosting after the unit has been turned off.

Serial # : _____

Install date : ____ / ____ / ____

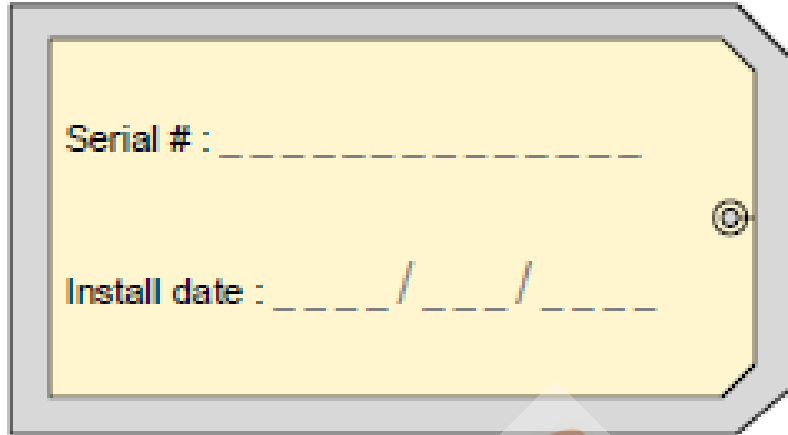

21750 Coolidge Highway • Oak Park, MI 48237 USA
+1 800-362-8491 • 248-291-2040 • Email: info@mopec.com

TABLE OF CONTENTS

UNPACKAGING YOUR PRODUCT.....4

INSTALLATION.....5

INTRODUCTION5

PRODUCT DRAWING

MB1006

SPECIFICATIONS.....7

OPERATING INSTRUCTIONS.....7

CLEANING AND MAINTENANCE

Stainless Steel Care and Maintenance.....7

Parts.....11

WARRANTY.....12

SAFETY LABELS.....13

TROUBLE SHOOTING GUIDE.....14

ELECTRICAL DIAGRAM.....15




UNPACKING

- 1) Carefully inspect the exterior of the shipping container before opening. If the crate is damaged and the product has sustained damage then immediately contact Mopec and the freight carrier. **Never discard the shipping container even if it is damaged beyond recognition.**
- 2) Have the delivery driver note any suspected damage on the Bill of Lading and sign it. Mopec will help assist in filing a claim for product repair and/or replacement.
- 3) Carefully open the containers and inspect the equipment for concealed damage. If visible damage is noticed (i.e. broken welds, dented stainless, scratches, etc.) follow through as noted above. Do not discard the shipping material. They are important in settling claims.

CAUTION: There are loose components in the packaging of your product. Be very careful in examining the packaging material as it may contain installation parts and/or product components.

INSPECTION

- 1) After carefully unpacking your Mopec Grossing station please inspect the items in the list below prior to installing the unit. The Grossing station should be thoroughly checked for loose screws, defects, or damage that may have occurred during shipping or packaging.

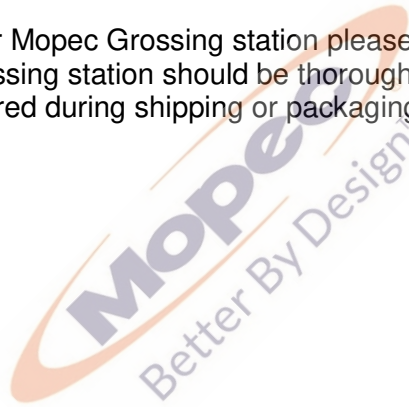

INSTALLATION

The BK105 chill tray is designed to work in your standard laboratory work area. The only requirement is, the unit must have a **minimum of 12" clearance to the front and back** louver panels. This allows the compressor to receive proper airflow for the compressor and exhaust.

If the BK105 Chill Tray was stored, or was laid in any position other than right side up, the unit must sit for 24 hours to allow the oil in the compressor to return to the operating position **BEFORE** plugging in and operating the unit.




INTRODUCTION

Thank you for purchasing the Mopec BK105 Chill Tray. With proper care and treatment, it will provide years of reliable service. It has been carefully checked at the factory and the tray pan temperature is pre-set at -20°F , + or - 3 degrees C

UNIT DRAWINGS

MB100

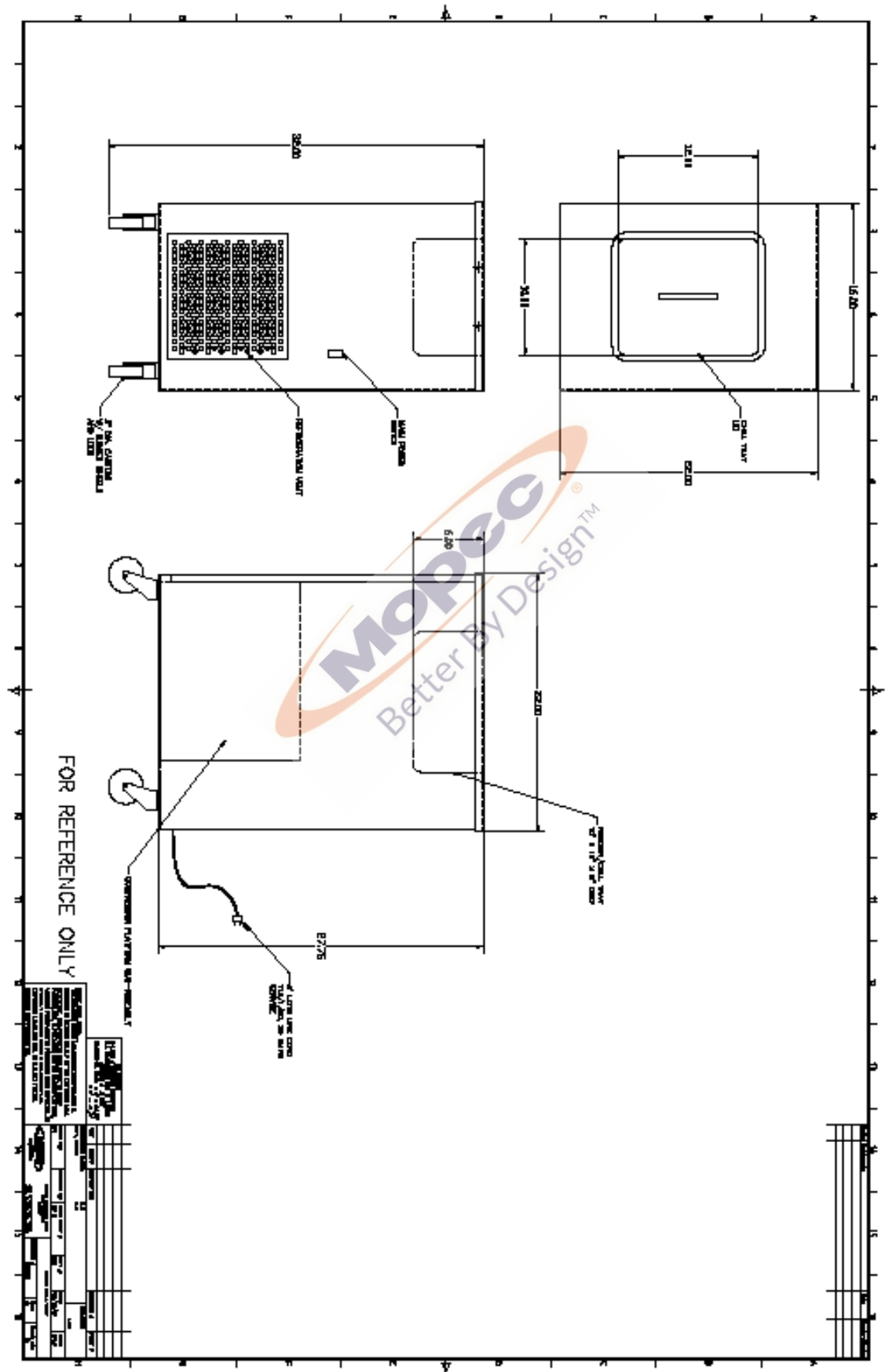

SPECIFICATIONS

- All Stainless-Steel Construction
- Casters: 3" Diameter and Swivel, Front 2 (two) Casters Locking
- Storage Chamber: 9 ¼" x 11 ½" x 6" deep. Complete with Cover
- Outside Dimensions: 16 ½" wide x 22 ½" deep x 32" high
- On/Off Rocker Switch
- 8 Foot Line Cord with Hospital Grade Plug

Electrical:	115 volts	60Hz	6 amps
Weight:	89 Pounds (net)		
Refrigeration:	R404A	Environmentally Safe	

OPERATION

1. Plug into 115-volt receptacle (8-foot line cord supplied on unit).
2. Push rocker switch on side of unit up to the on position. The compressor will start. Tray temperature factory setting is at -20 degrees +/- 3 degree. Please allow 30 to 45 minutes for unit to reach temp. If you need to adjust the temperature of the unit, then please call Mopec Service at (800) 362-8491. Our technician will walk you through the proper steps to make this adjustment.

AIR HANDLING

The front and back of the BK105 must be kept a minimum of 12 inches from a obstruction. This will provide enough air flow to keep the nit operating efficiently and without icing up from lack of air circulation.

CLEANING AND MAINTENANCE

DISINFECTING STAINLESS STEEL SURFACES

All stainless steel surfaces can be cleaned with soap and water, which will remove debris.

The stainless steel surfaces can be disinfected with a non-caustic disinfectant. We suggest using BE045 Path Cloud or BE047 Bench Wipe for cleaning purposes. We recommend you **NOT USE** a bleach solution to clean your unit. Bleach will eventually erode stainless steel if not thoroughly rinsed. **The use of chlorine bleach will VOID THE STAINLESS STEEL WARRANTY**

To maintain your stainless-steel product, follow these four steps:

1) **Never, ever use wire brushes, Brillo, steel wool or abrasive cleansers (like Ajax or Comet).** If something needs to be aggressively cleaned only use a Scotch-brite pad or similar product and only scour with the "grain" of the stainless. As an example, please reference the photo on page 2 of this document. It is clear that a very abrasive product was used in an area on the unit and that did not follow the grain of the stainless. The effects of this scratching may diminish over time with proper care but the effects of this scouring are

obvious. (The use of the Scotch-brite Pad following the grain over time may help scratches such as this). Depending on the surface finish of your stainless steel, abrasive cleaners can cause scratching. Duller finishes probably won't show scratching as much as mirror or highly polished finishes. When in doubt, test in a hidden spot, and also work from the least risky type of cleaning, (i.e. water) to the heavy-duty stuff.

Do not use cleaners containing chlorine. While it may be second nature to bleach everything, stainless steel and chlorine do not mix well. Do not use bleach when cleaning stainless steel. Do not allow bleach or bleach water to sit for long periods. Bleach can eventually cause staining and pitting. Bleach stains are removed with stainless steel cleaning polish.

2) Keep the surface clean of grime, tissue and particulates. This can be accomplished by using the "Water Hand Spray Rinse" and use of cleaning products.

3) Rinse the surface after using disinfectant. In Pathology and other medical areas, the act of disinfecting is desired. There are a number of ways to do this including using Mopec's Bench Spray & Wipe Disinfectant. Labs use any number of different products including 10% bleach or other disinfectant sprays and wipes. For the most part, each and every one of these has high salt contents and lower PH levels to aid with disinfection. Most disinfectants must be followed up with a water rinse to remove the salts that remain after these products dry. We advise to always follow up a disinfection cleaning with a thorough rinse of water. We advise not to use diluted bleach, if you must; we stress the importance of a thorough rinse of water after use. If not rinsed properly, these salts can become visible after the disinfectant dries. They can appear with a whitish characteristic or contain light lines of white with a grainy feel when you wipe your hand across the work surface. If these residues are not removed with a thorough water rinse and wipe down they will accumulate and eventually degrade the appearance and integrity of your stainless surface. Rust is a long-term possibility if salts are allowed to remain on the work surfaces over time. Gritty, dirty water or residue from cleaning solutions left on a stainless-steel surface can stain or damage the finish.

4) All stainless-steel products should be protected by a polish. As a prime example before any product leaves Mopec it is coated with WD-40 as a protecting coating for the stainless. Mopec offers a Stainless-Steel Cleaner and Polish in both wipe and spray. These Mopec products will not only deep clean your stainless but will also protect their finish from chemical, low PH and salt and keep the finish looking like new.

Decal

Mopec advises that if the technicians are not taking the proper precautions when using the Decal solution, the possibility of two things will occur, a brown or rust ring where the Decal resides along with a milky white substance on the surface. (As an example, reference the below photos to see the rust rings.) Decal is very harsh, even the fumes can and will cause staining on stainless steel. One thing that you might want to consider doing is to place the Decal container you currently use inside a plastic base that will help catch drips that might occur. Clean and rinse your station after every use of Decal solution.

Rust

Rust can and will occur on stainless if it is not maintained properly. The most common cause of rust is from using metal or stainless racks that are not made of 304 stainless. This is referred to as "transfer rust". Leaving of salts from cleaners or disinfectants can and will lead to possible rust areas in the long term. Always rinse all disinfectants before they dry. Decal solutions and even fumes are very aggressive and can cause rust if not cleaned up and used properly around stainless. Formalin use has not been shown to cause rust in any way.

Conclusions & Suggestions

We are confident and can assure you that if you institute the suggestions detailed above that your Mopec Grossing station will look as it did the day it arrived.



Do not assume it's the cleaner. If you do have some spotting or staining, and you've followed all of the suggestions, it may not be the cleaner. Water, especially hard water, can leave spotting and staining on stainless steel surfaces. Hard water can leave mineral deposits, resulting in whitish-colored spots and streaks. Remove hard water stains with vinegar or with stainless steel cleaning polish. Prevent hard water stains by towel-drying after every wash.

Do not allow soaps and cleaners to dry on surfaces. The chemicals in many soaps and cleaners can cause staining. Never use corrosive cleaners such as mineral spirits. Use stainless steel cleaning polish and a

non-abrasive scrub pad to remove dried cleaner stains. Baking soda mixed with liquid dish soap can make a good paste to gently rub on stains. Be sure to rinse the stainless-steel surface thoroughly, and towel dry.

If stains remain Mopec recommends trying a stainless-steel cleaner and polisher. Barkeeper's Friend is a good powder formula that can clean without scratching. Be sure to follow the directions, rinse thoroughly, and towel dry. These methods should help remove any discoloration.

Fingerprints and Stains – The most common surface contaminants that occur from normal use are fingerprints and mild stains. These usually affect only appearance so fortunately they do not have an effect on corrosion resistance. They can easily be removed by a variety of simple cleaning methods. The most troublesome marks to remove from the surface of smooth polished or bright finished stainless steel are fingerprints; fortunately, they can be removed with a common glass cleaner or by gently rubbing with a paste of soda ash (sodium carbonate) and water which would be applied with a soft cloth. Again, it is best to follow with a warm water rinse.

Clean Water and Wipe – The method that will do an adequate job and is the simplest, safest and the least costly is the best method. There is no surface coating to wear off of stainless steels so the surface will thrive with frequent cleaning. The first choice to clean mild stains and loose dirt and soil should always be a soft cloth and clean, warm water. Rinsing with clean water and wiping the surface dry will finish the process and eliminate the possibility of water stains.

Solvent Cleaning – To remove oils, greases and fresh fingerprints that have not had time to oxidize or decompose, use a solvent that does not contain chlorine. There are many organic cleaners on the market today that optimize safety attributes and clean ability. Spray or vapor methods or by wiping with clothes containing solvents can also clean surfaces. The wiping technique will sometimes leave the surface streaked.

Scratch Repair

A surface scratch can be repaired using the following technique. Completely removing the scratch will depend on how severe it is.

Use 120 grit emery cloth or paper and firm pressure to sand the scratch. Sanding must always go in the direction of the grain. Sand in a perfectly straight line, avoiding the natural tendency to sand in an arc. Sand the surface until the scratch is gone.

Polish using a very fine grade of 3M scotch-brite pads. Use the same motions as with sanding. Polish the surface until the original finish is restored.

For stubborn spots, stains, light discoloration, water marking or light rust staining use a mild, non-scratching cream or polish. Apply with soft cloth or soft sponge and rinse off residues with clean water and dry. Avoid cleaning pastes with abrasive additions. Suitable cream cleansers are available with soft calcium carbonate additions, or with the addition of citric acid. Do not use chloride or acidic solutions



Nylon abrasive pads should be adequate for dealing with most deposits (DO NOT USE STEEL WOOL OR BRILLO PADS). If a more severe treatment is needed to mask coarse scratches or physical damage on a surface, use the finest abrasive medium consistent with covering the damage marks. With directional brushed and polished finishes, align and blend the new "scratch pattern" with the original finish, checking that the resulting finish is aesthetically acceptable. Silicon carbide media may be used, especially for the final stages of finishing. Avoid using hard objects such as knife blades and certain abrasive/souring agents as it is possible to introduce surface scuffs and scratches. Scratching is particularly noticeable on sink drainer areas. These are usually superficial and can be removed with proprietary stainless steel cleaners or, alternatively, with a car paint restorer, such as 'T-cut'. Rust marks or staining on stainless steels is unlikely to be the result of corrosion to the stainless steel itself (similar marks may also be found on porcelain and plastic sinks). These marks are likely to result from small particles of carbon steel from wire wool.

Replacement parts are available from Mopec The can be ordered by contacting Mopec at 800-362-8491.

BK105 Chill Tray Parts List

<i>Blank Cover - Handy Box</i>	<i>2MULBLC-CW</i>
<i>Caster</i>	<i>PD0002</i>
<i>Caster W/Lock</i>	<i>PD0003</i>
<i>Compressor</i>	<i>2M4FCOM-CW</i>
<i>Cover, Pan Steam</i>	<i>PM0349</i>
<i>Defrost Timer</i>	<i>EDT10</i>
<i>Drier</i>	<i>2C03DRI-CW</i>
<i>Front Grill</i>	<i>PF1050</i>
<i>Handy Box</i>	<i>2EGSHAB-CW</i>
<i>Pan, Stainless Steel</i>	<i>PM0346</i>
<i>Pipe Insulation</i>	<i>2612INS-CW</i>
<i>Pipe Insulation</i>	<i>2I38INS-CW</i>
<i>R404 Tag</i>	<i>2404TAG-CW</i>
<i>Refrigerant 404A</i>	<i>2404RIF-CW</i>
<i>Switch w Pilot Light</i>	<i>PE0037</i>
<i>Temperature Control</i>	<i>060H1101 Danfoss</i>

LIMITED WARRANTY

Products manufactured by Mopec will be free from defects in material and workmanship and conform to Mopec's description or specifications. If a warranty claim is made within one (1) year from the earlier of the date of acceptance/first beneficial use, the defective or nonconforming Product or Part thereof will be repaired or (at Mopec's option) replaced free of charge, FCA Mopec's dock. All warranty claims must be in writing and received by Mopec within the warranty period. The warranty is not transferable (other than to customers of Mopec's authorized Distributors), and will not apply unless the Equipment has been properly installed, maintained and operated in accordance with all instructions; and does not apply to defects, nonconformities or other failure due to Equipment misuse, abuse, modifications, or other causes outside Mopec's control. If a warranty claim is made in writing within the warranty period, the defective or nonconforming Equipment (or Part thereof) will be repaired or (at Mopec's option) replaced free of charge, FCA Mopec's dock.


THERE ARE NO WARRANTIES WHICH EXTEND BEYOND THE DESCRIPTION ON THE FACE HEREOF. THE WARRANTY AS SET FORTH HEREIN IS IN LIEU OF ALL OTHER WARRANTIES, EXPRESS OR IMPLIED, INCLUDING ANY IMPLIED WARRANTY OF MERCHANTABILITY OR FITNESS FOR A PARTICULAR PURPOSE.

To the extent that Mopec is acting as a supplier of Products manufactured by a third party, the Products will be warranted only to the extent that they are warranted by their manufacturers and Buyer (or its customer) agrees to look solely to the Product manufacturer for all warranty claims.

MOPEC, 21750 COOLIDGE HIGHWAY, OAK PARK, MI 48237

SAFETY LABELS

BEFORE SERVICING THE UNIT LOOK FOR AND HEED THE FOLLOWING LABEL

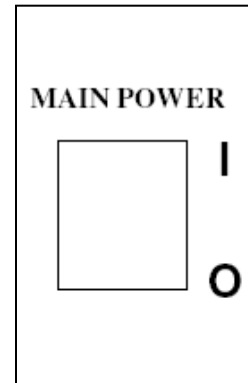

TROUBLE SHOOTING

Power Switch is on and nothing happens

1. Check the Unit to ensure it is plugged in
2. Check the circuit breaker for the outlet to ensure the breaker is not tripped
3. If your BK105 Chill Tray does not come on when you first turn the switch, leave the unit on and plugged in for approximately 45 minutes. The unit may have entered its defrost cycle when the unit was shut off. The defrost timer will not let it start the cooling cycle until the timer has completed the cycle. Once the defrost cycle has been completed the unit will start to cool down.
4. If the unit has been plugged in for more than an hour and still has not started to chill call customer service at 800-362-8491


The Unit does not reach the required temperature

1. The room conditions may be too humid or too hot for the optimum operating conditions. High heat and or High Humidity will adversely affect the unit's ability to cool.
2. The evaporator coil may need to be cleaned to allow proper air flow.

Water on floor after turning the unit off

1. It is normal for small amounts of water to build up underneath the unit due to condenser defrosting.

ELECTRICAL DIAGRAM

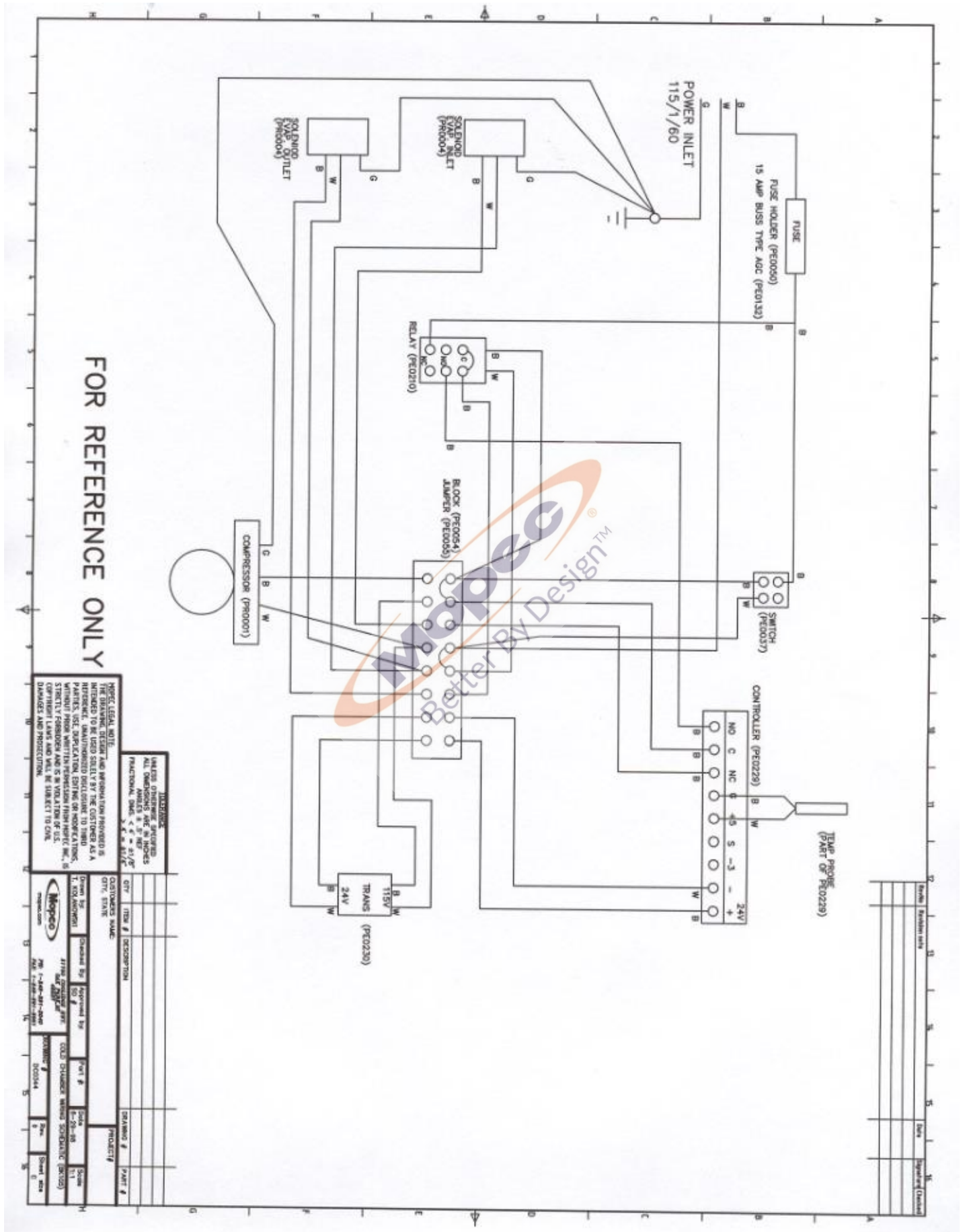

FOR REFERENCE ONLY

WARRANTY INFORMATION

THESE TERMS AND CONDITIONS APPLY TO ALL MOPCO PRODUCTS. ALL DIMENSIONS ARE IN INCHES UNLESS OTHERWISE SPECIFIED. MOPCO SHALL NOT BE LIABLE FOR DAMAGES AND LOSS OF PROFITS.

WARRANTY: MOPCO warrants its products to be free from defects in material and workmanship for a period of 12 months from the date of purchase. This warranty is void if the product has been used in an unauthorized manner or if the product has been damaged by fire, flood, or other causes beyond Mopco's control.

REPAIRS: Repairs shall be made at Mopco's expense if the product is found to be defective within the warranty period. Repairs shall be made at the customer's expense if the product is found to be defective after the warranty period or if the product has been damaged by fire, flood, or other causes beyond Mopco's control.

CLAIMS: Claims shall be made in writing to Mopco within 30 days of the date of purchase. Mopco shall not be liable for claims for consequential or incidental damages.

FORCE MAJEURE: Mopco shall not be liable for delays or non-performance due to acts of God, war, strikes, or other causes beyond Mopco's control.

ASSIGNMENT: These terms and conditions shall apply to all assignments of the product.

ENTIRE AGREEMENT: These terms and conditions constitute the entire agreement between Mopco and the customer.

GOVERNING LAW: These terms and conditions shall be governed by the laws of the state of California.

CONTACT: Mopco, 11111 S. Bascom Ave., Suite 100, San Jose, CA 95128. Phone: (408) 261-1111. Fax: (408) 261-1112. Website: www.mopco.com



Better By Design™

21750 Coolidge Highway • Oak Park, MI 48237 USA
+1 800-362-8491 • 248-291-2040 • Email: info@mopec.com

2015 Mopec all rights reserved

