




USER MANUAL

Poly Body Slats BC011



Serial # : _____

Install date : ____ / ____ / ____

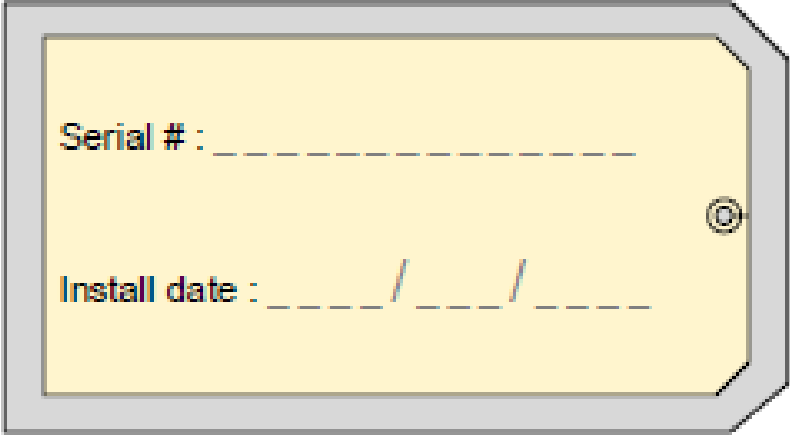

21750 Coolidge Highway • Oak Park, MI 48237 USA
+1 800-362-8491 • 248-291-2040 • Email: info@mopec.com

TABLE OF CONTENTS

UNPACKAGING YOUR PRODUCT.....4

INSTALLATION.....4

INTRODUCTION4

CLEANING AND MAINTENANCE

Disinfecting and Cleaning.....5

WARRANTY.....6

UNPACKING

- 1) Carefully inspect the exterior of the shipping container before opening. If the crate is damaged and the product has sustained damage then immediately contact Mopec and the freight carrier. **Never discard the shipping container even if it is damaged beyond recognition.**
- 2) Have the delivery driver note any suspected damage on the Bill of Lading and sign it. Mopec will help assist in filing a claim for product repair and/or replacement.
- 3) Carefully open the containers and inspect the equipment for concealed damage. If visible damage is noticed (i.e. broken welds, dented stainless, scratches, etc.) follow through as noted above. Do not discard the shipping material. They are important in settling claims.

CAUTION: There are loose components in the packaging of your product. Be very careful in examining the packaging material as it may contain installation parts and/or product components.

INSPECTION

- 1) After carefully unpacking your Mopec Body Slats please inspect the items in the list below prior to installing the unit.

INSTALLATION

To install the BC011 body slats simply place under the body in the desired locations to elevate the body of the tray or surface to allow better draining of fluids. The long flat side is placed against the body: typical placement are shoulders, small of back, buttocks and legs.

INTRODUCTION

Set of four body slats elevate the body on a table during the autopsy or mortuary preparation. While elevated, waste drains without obstruction. Suggested placement of slats are head, shoulders, buttocks and heels. Made of a chemically resistant white polyethylene, each measuring almost 3" wide (7.6 cm) and over 22" long (57.5 cm), can be cleaned using water and one of Mopec's specially formulated cleaners.

Designed for use in the 23" JC100 Body Tray but can be used separately.

CLEANING AND MAINTENANCE

DISINFECTING AND CLEANING STAINLESS STEEL SURFACES

All stainless steel surfaces can be cleaned with soap and water, which will remove debris.

The stainless steel surfaces can be disinfected with a non-caustic disinfectant. We suggest using BE045 Path Cloud or BE047 Bench Wipe for cleaning purposes. We recommend you **NOT USE** a bleach solution to clean your unit. Bleach will eventually erode stainless steel if not thoroughly rinsed. **The use of chlorine bleach will VOID THE STAINLESS STEEL WARRANTY**

During the cleaning we suggest wiping the surface in the same direction as the satin finish which will help lift up dirt from the grain finish.

Most scratches can be removed simply by utilizing a "non-metallic" abrasive pad and rubbing in the same direction as the satin finish.

Since most abrasive pads vary from one supply to another, we suggest rubbing the entire surface to blend the scratch and blend the balance of the surface.

Plexi-glass surfaces will scratch if cleaned with an unsuitable cleaner and improper cleaning. Many plastic cleaners are available and we suggest using one. Wipe dry with a clean, absorbent cloth or paper towel turning often

LIMITED WARRANTY

Products manufactured by Mopec will be free from defects in material and workmanship and conform to Mopec's description or specifications. If a warranty claim is made within one (1) year from the earlier of the date of acceptance/first beneficial use, the defective or nonconforming Product or Part thereof will be repaired or (at Mopec's option) replaced free of charge, FCA Mopec's dock. All warranty claims must be in writing and received by Mopec within the warranty period. The warranty is not transferable (other than to customers of Mopec's authorized Distributors), and will not apply unless the Equipment has been properly installed, maintained and operated in accordance with all instructions; and does not apply to defects, nonconformities or other failure due to Equipment misuse, abuse, modifications, or other causes outside Mopec's control. If a warranty claim is made in writing within the warranty period, the defective or nonconforming Equipment (or Part thereof) will be repaired or (at Mopec's option) replaced free of charge, FCA Mopec's dock.

THERE ARE NO WARRANTIES WHICH EXTEND BEYOND THE DESCRIPTION ON THE FACE HEREOF. THE WARRANTY AS SET FORTH HEREIN IS IN LIEU OF ALL OTHER WARRANTIES, EXPRESS OR IMPLIED, INCLUDING ANY IMPLIED WARRANTY OF MERCHANTABILITY OR FITNESS FOR A PARTICULAR PURPOSE.

To the extent that Mopec is acting as a supplier of Products manufactured by a third party, the Products will be warranted only to the extent that they are warranted by their manufacturers and Buyer (or its customer) agrees to look solely to the Product manufacturer for all warranty claims.

MOPEC, 21750 COOLIDGE HIGHWAY, OAK PARK, MI 48237



21750 Coolidge Highway • Oak Park, MI 48237 USA
+1 800-362-8491 • 248-291-2040 • Email: info@mopec.com

2015 Mopec all rights reserved