

HC Series

HC100 HC150 HC200
HC250 HC300 HC350 HC375
HOODED DISSECTING TABLES



REFERENCE MANUAL



21750 Coolidge Highway • Oak Park, MI 48237 USA
+1 800-362-8491 • 248-291-2040 • Email: info@mopec.com

IMPORTANT INFORMATION


MODEL NO.:

SERIAL NO.:

DATE OF PURCHASE:

DATE OF INSTALLATION:

INSTALLATION NOTES:

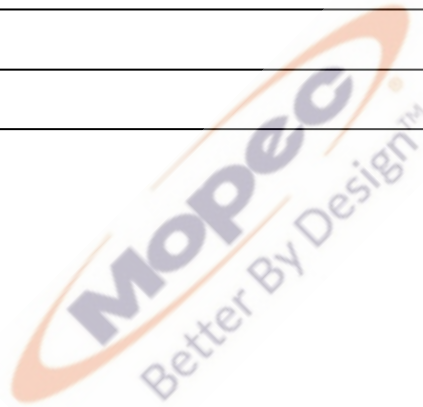

SERVICE AND SUPPORT:



21750 Coolidge Highway • Oak Park, MI 48237 USA
+1 800-362-8491 • 248-291-2040 • Email: info@mopec.com

TABLE OF CONTENTS

INTRODUCTION	4-6
UNPACKAGING YOUR PRODUCT	7
CLEANING AND MAINTENANCE	8
WARRANTY	9
PRODUCT DRAWINGS	10

HC100 Rough In	10
HC100	11
HC150 Rough In	12
HC150	13
HC200 Rough In	14
HC200	15
HC250 Rough In	16
HC250	17
HC300 Rough In	18
HC300	19
HC350 Rough In	20
HC350	21
HC375 Rough In	22
HC375	23

INTRODUCTION

The Mopec HC Series Dissection Tables with ventilation hood is a versatile and economical dissecting work system for pathology, histology and necropsy laboratories.

The table is designed to be connected to existing ventilation systems to provide a efficient and maintenance-free ventilated workstation.

HC100 Dissection Table with Vented Hood - 60" Right Sink

RIGHT SINK

- All stainless steel construction
- Dimensions: 60" Long x 30" Wide x 37" High Work area
- Overall unit height, 55"
- Two (2) low profile fluorescent light fixtures
- Full length Lexan splash shield
- Large deep sink located on right
- Hot/Cold water fixture with wrist blade handles & swing spout faucet (vacuum breaker protected)
- GFCI Electrical receptacle
- Vent tie-in, 36" L x 4" W
- All internal plumbing and electrical lines are factory installed and require only trade hook-up

HC150 Dissection Table with Vented Hood - 60" Left Sink

LEFT SINK

- All Stainless Steel Construction
- Dimensions: 60" L x 30" W x 37" H work area.
- Overall unit height, 55"
- Two (2) low profile fluorescent light fixtures
- Full length Lexan splash shield
- Large deep sink located on left
- Hot/Cold water fixture with wrist blade handles & swing spout
- GFCI Electrical receptacle
- Vent tie-in 36" L x 4" W
- All internal plumbing and electrical lines are factory installed and require only trade hook-up

HC200 Dissection Table with Vented Hood - 72" Right Sink

RIGHT SINK

- All stainless steel construction
- Dimensions: 72" Long x 30" Wide x 37" High Work Area
- Overall Height, 55"
- Two (2) low profile fluorescent light fixtures
- Full length Lexan splash shield
- Large deep sink located on right
- Hot/Cold water fixture with wrist blade handles & swing spout faucet (vacuum breaker protected)
- GFCI Electrical Receptacle
- Vent tie-in, 48" L x 4" W
- All internal plumbing and electrical lines are factory installed and require only trade hook-up

HC250 Dissection Table with Vented Hood - 72" Left Sink

LEFT SINK

- All Stainless Steel Construction
- Dimensions: 72" Long x 30" Wide x 37" High work area
- Overall unit height, 55"
- Two (2) low profile fluorescent light fixtures
- Full length Lexan splash shield
- Large deep sink located on left
- Hot/Cold water fixture with wrist blade handles & swing spout faucet (vacuum breaker protected)
- G.F.C.I. Electrical receptacle
- Vent tie-in, 48" L x 4" W
- All internal plumbing and electrical lines are factory installed and require only trade hook-up

HC300 Dissection Table with Vented Hood - 84" Right Sink

RIGHT SINK

- All stainless steel construction
- Dimensions: 84" Long x 30" Wide x 37" High work area
- Overall height, 55")
- Two (2) low profile fluorescent light fixtures
- Full length Lexan splash shield
- Large deep sink located on right
- Hot/Cold water fixture with wrist blade handles & swing spout faucet (vacuum breaker protected)
- G.F.C.I. Electrical receptacle
- Vent tie-in 60" L x 4" W
- All internal plumbing and electrical lines are factory installed and require only trade hook-up

HC350 Dissection Table with Vented Hood - 84" Left Sink

LEFT SINK

- All stainless steel construction
- Dimensions: 84" Long x 30" Wide x 37" High work area
- Overall height, 55"
- Two (2) low profile fluorescent light fixtures
- Full Length Lexan Splash Shield
- Large deep sink located on left
- Hot/Cold water fixture with wrist blade handles & swing spout faucet (vacuum breaker protected)
- G.F.C.I. Electrical receptacle
- Vent tie-in, 60" L x 4" W
- All internal plumbing and electrical lines are factory installed and require only trade hook-up

HC375 Dissection Table with Vented Hood - 84" Center Sink

CENTER SINK

- All Stainless Steel Construction
- Dimensions: 84" Long x 30" Wide x 37" High work area
- Overall height, 55"
- Two (2) low profile fluorescent light fixtures
- Full length Lexan splash shield
- Large deep sink located in the center of unit
- Hot/Cold water fixture with wrist blade handles & swing spout faucet (vacuum breaker protected)
- G.F.C.I. Electrical receptacle
- Vent tie-in, 60" L x 4" W
- All internal plumbing and electrical lines are factory installed and require only trade hook-up

Available Options

HO007: Hand Spray Rinse

HO008: Gas Valve Deck Mounted

HO009: Air Valve Deck Mounted

BL800: Disposal 1 H.P. Heavy Duty

UNCRATING / SET-UP

The first step in using your new Dissecting Table is to carefully inspect the exterior of the shipping container before opening. If the crate is damaged and the product has sustained damage than immediately contact Mopec and the carrier and never discard the shipping container even if they're damaged beyond recognition.

The receiver must note any suspected damage on the bill of lading and sign it. Mopec will help assist in filing a claim for product repair and/or replacement.

Carefully open the containers and inspect the goods for concealed damage. If concealed damage is noticed (i.e. cracked Plexiglas, broken welds, dented stainless, etc.) follow through as noted above being certain not to discard packing materials.



CLEANING INSTRUCTIONS

DISINFECTING AND MAINTAINING STAINLESS STEEL SURFACES

All stainless steel surfaces can be cleaned with soap and water, which will remove debris.

The stainless steel surfaces can be disinfected with a non caustic disinfectant. We suggest using BE045 PathCloud or BE047 Bench Wipe for cleaning purposes.

We recommend you **DO NOT USE** a bleach solution to clean your unit, bleach will eventually erode stainless steel if not thoroughly rinsed. **The use of chlorine bleach will VOID THE STAINLESS STEEL WARRANTY**

During the cleaning we suggest wiping the surface in the same direction as the satin finish which will help elevate dirt from the grain finish.

Scratches can be removed simply by utilizing a "non-metallic" abrasive pad and rubbing in the same direction as the satin finish.

Since most abrasive pads vary from one supply to another, we suggest rubbing the entire surface to blend the scratch and blend the balance of the surface.

MOPEC DISSECTING TABLE WARRANTY

DISSECTING TABLE EQUIPMENT AND ALL PARTS THEREOF, MANUFACTURED BY MOPEC IS WARRANTED TO OUR CUSTOMERS TO BE FREE FROM DEFECTS IN MATERIAL AND CONSTRUCTION WHEN SUBJECTED TO NORMAL USE AND SERVICE. THIS WARRANTY SHALL NOT APPLY TO ANY EQUIPMENT OR PART THEREOF, WHICH HAS BEEN SUBJECT TO ALTERATIONS, ACCIDENT, MISUSE, OR WHICH HAS BEEN USED MORE THAN ITS PUBLISHED CAPACITY.

OUR OBLIGATION UNDER THIS WARRANTY IS LIMITED TO REPAIRING OR REPLACING, WITHOUT CHARGE ANY PART OR PARTS OF SAID EQUIPMENT WHICH PROVE DEFECTIVE AND WHICH OUR EXAMINATION SHALL DISCLOSE TO OUR SATISFACTION TO BE THUS DEFECTIVE. WE DO NOT ACCEPT RESPONSIBILITY FOR CONSEQUENTIAL DAMAGE RESULTING FROM SUCH DEFECTIVE PARTS. THE PERIOD OF WARRANTY ON MOPEC EQUIPMENT IS AS FOLLOWS:

STAINLESS STEEL MATERIAL: STAINLESS STEEL IS WARRANTED FOR LIFE, TO BE FREE OF CORROSION. THIS WARRANTY IS APPLICABLE ONLY IF THE MATERIAL IS CLEANED AND DISINFECTED WITH A CLEANER THAT IS NOT CAUSTIC AND APPROVED FOR STAINLESS. SEE CLEANING AND DISINFECTING PROCEDURES.

ALL EQUIPMENT: EQUIPMENT IS WARRANTED IN ITS ENTIRETY FOR A PERIOD OF ONE (1) YEAR FROM THE DATE OF SALE. THIS WARRANTY INCLUDES PARTS AND LABOR FOR COVERED REPAIRS.

MOPEC WILL ACCEPT INCOMING DEFECTIVE PARTS, ONLY WHEN SHIPMENT IS PREPAID. THIS WARRANTY CONTRACT IS IN LIEU OF ALL OTHER WARRANTIES, AND RELEASES MOPEC OF ALL OTHER OBLIGATIONS OR LIABILITIES. IT NEITHER ASSUMES NOR AUTHORIZES ANY PERSON OR PERSONS TO ASSUME ANY OBLIGATION OTHER THAN THAT COVERED IN THIS WARRANTY.

MOPEC 21750 COOLIDGE HIGHWAY OAK PARK, MI 48237

