

OA300

NECROPSY/DISSECTING TABLE



REFERENCE MANUAL



21750 Coolidge Highway • Oak Park, MI 48237 USA
+1 800-362-8491 • 248-291-2040 • Email: info@mopec.com

IMPORTANT INFORMATION

MODEL NO.:


SERIAL NO.:

DATE OF PURCHASE:

DATE OF INSTALLATION:

INSTALLATION NOTES:

SERVICE AND SUPPORT:

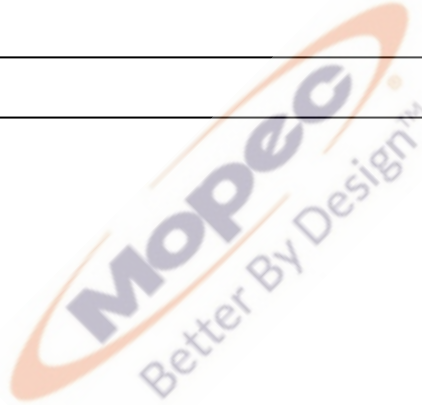

21750 Coolidge Highway • Oak Park, MI 48237 USA
+1 800-362-8491 • 248-291-2040 • Email: info@mopec.com

TABLE OF CONTENTS

INTRODUCTION	Page 4
UNCRATING	Page 5
INSTALLATION	Page 5
CLEANING INSTRUCTIONS	Page 6
WARRANTY	Page 7



INTRODUCTION

The Mopec OA300 Necropsy/Dissection Table is an economical workstation ideal for small and large laboratories. Standard design features include the sloped work area, interchangeable grid plates. Below is a brief list of its features:

- The tabletop is fabricated of 14 Ga type stainless steel finished with a #4 satin finish, with integral sinks.
- Swing spout faucets are provided with vacuum breaker and wrist blade H/C handles, commercial quality cast brass chrome plated.
- Hand Held Table Rinse / Spray which is deck mounted for convenience and are fabricated of brass chrome plated along with high pressure nylon reinforced tubing.
- Perforated Grid Plates: Fabricated of 16 Gauge each plate is perforated with 1/2" perforations on 1" centers, each grid panel is self supporting with 1/2" recess and completely interchangeable. (2) Grid plates, all interchangeable.
- All internal Plumbing lines are factory installed and require only trade hook-up.

Available Option:

- Heavy-Duty Heavy Duty Disposal 1 HP, 115V/60Hz. Complete with on/off switch, water solenoid and vacuum breaker plumbed and wired ready for use.

UNCRATING

The first step in installing your new Necropsy Table is to carefully inspect the exterior of the shipping container before opening. If the crate is damaged and the product has sustained damage than immediately contact Mopec and the carrier, never discard the shipping container even if they're beyond recognition. Have the driver note any suspected damage on the bill of lading and sign it. Mopec will help assist in filing a claim for product repair and/or replacement. Carefully open the containers and inspect the goods for concealed damage. If concealed damage is noticed (i.e. cracked Plexiglas, broken welds, dented stainless, etc.) following through as noted above, and again does not discard the shipping material, they are important in settling claims.

INSTALLATION

Once the necropsy sink is unpacked, carefully place it in the area where it is to be used. Align the edge of the table such that it's true and positioned according to the facility requirements. Level the table by placing a level on the table edges both sides to side and end to end do not place the level on the table surface. "Important", the table surface is sloped toward the drain and will yield incorrect results.

Once the sink is leveled, the utilities can be hooked up. The cold water line is 1/2" and the hot water is 1/2" hard copper which will require a customer supplied on/off valves. We suggest ball type water valves to assure proper water flow in operating all table accessories. Drain lines are standard 1-1/2" IPS and need to be vented to assure proper drainage. Electrical requirements are solely dependent on accessories. If the table excludes a disposal than one 115/1/60, 15 amps circuit is sufficient. If a disposal is included that another 115/1/60, 15 amp circuits must be added, this is tied in to the same junction box yet wired independently.

CLEANING INSTRUCTIONS

All stainless steel surfaces can be cleaned with soap and water, which will remove debris. The stainless steel surfaces can be disinfected with a not caustic disinfectant. A 10% bleach solution can be used but must be rinsed very well; bleach will eventually erode stainless steel if not rinsed. During the cleaning we suggest wiping the surface in the same direction as the satin finish which will help elevate dirt from the grain finish. Scratches can be removed simply by utilizing a "non-metallic" abrasive pad and rubbing in the same direction of the satin finish. Since most abrasive pads vary from one supply to another, so we suggest rubbing the entire surface to blend the scratch and blend the balance of the surface.



MOPEC NECROPSY TABLE WARRANTY

NECROPSY TABLE EQUIPMENT AND ALL PARTS THEREOF, MANUFACTURED BY MOPEC, AND IS WARRANTED TO OUR CUSTOMERS TO BE FREE FROM DEFECTS IN MATERIAL AND CONSTRUCTION WHEN SUBJECTED TO NORMAL USE AND SERVICE. THIS WARRANTY SHALL NOT APPLY TO ANY EQUIPMENT OR PART THEREOF, WHICH HAS BEEN SUBJECT TO ALTERATIONS, ACCIDENT, MISUSE, OR WHICH HAS BEEN USED MORE THAN ITS PUBLISHED CAPACITY.

OUR OBLIGATION UNDER THIS WARRANTY IS LIMITED TO REPAIRING OR REPLACING, WITHOUT CHARGE ANY PART OR PARTS OF SAID EQUIPMENT WHICH PROVE DEFECTIVE AND WHICH OUR EXAMINATION SHALL DISCLOSE TO OUR SATISFACTION TO BE THUS DEFECTIVE. WE DO NOT ACCEPT RESPONSIBILITY FOR CONSEQUENTIAL DAMAGE RESULTING FROM SUCH DEFECTIVE PARTS. THE PERIOD OF WARRANTY ON MOPEC EQUIPMENT IS AS FOLLOWS:

STAINLESS STEEL MATERIAL: STAINLESS STEEL IS WARRANTED FOR LIFE, TO BE FREE OF CORROSION. THIS WARRANTY IS APPLICABLE ONLY IF THE MATERIAL IS CLEANED AND DISINFECTED WITH A CLEANER THAT'S NOT CAUSTIC AND APPROVED FOR STAINLESS. SEE CLEANING AND DISINFECTING PROCEDURES.

ALL EQUIPMENT: EQUIPMENT IS WARRANTED IN ITS ENTIRETY FOR A PERIOD OF ONE (1) YEAR FROM THE DATE OF SALE. THIS WARRANTY INCLUDES PARTS AND LABOR FOR COVERED REPAIRS.

MOPEC WILL ACCEPT INCOMING DEFECTIVE PARTS, ONLY WHEN SHIPMENT IS PREPAID. THIS WARRANTY CONTRACT IS IN LIEU OF ALL OTHER WARRANTIES, AND RELEASES MOPEC OF ALL OTHER OBLIGATIONS OR LIABILITIES. IT NEITHER ASSUMES NOR AUTHORIZES ANY PERSON OR PERSONS TO ASSUME ANY OBLIGATION OTHER THAN THAT COVERED IN THIS WARRANTY.

MOPEC, 21750 COOLIDGE HIGHWAY, OAK PARK, MI 48237.