

REV	DESCRIPTION	DATE	ECN #	APPR
A	INITIAL RELEASE	00-XXX-19	XXXXX	MM


THIS DOCUMENT CONTAINS PROPRIETARY INFORMATION OF MOPEC, AND MAY NOT BE COPIED IN WHOLE OR PART, AND NO USE SHALL BE MADE OF IT OTHER THAN FOR WHICH IT HAS BEEN SUBMITTED WITHOUT THE WRITTEN PERMISSION OF A MOPEC. MANAGER

MOPEC 2024

UNLESS OTHERWISE SPECIFIED:  
ALL DIMENSIONS IN INCHES

TOLERANCES	0.0	± 0.1 in
	0.00	± 0.05 in
	0.000	± 0.010 in
	0.0000	± 0.0010 in
	ANGLES	±0.5°


BREAK ALL CORNERS 0.063 X 45°  
ALL FILLET RADII R0.063 /0.125in  
INSPECT OVAL DIMENSIONS



MOPEC GROUP

subcontractor MANUFACTURING

DRAWN TO 3RD ANGLE PROJECTION

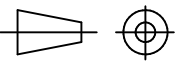

DRAWN BY: B. LAY  
DATE: 1/29/2020


CHECKED BY: \_\_\_\_\_  
DATE: 1/29/2020

APPROVED BY: \_\_\_\_\_  
DATE: M.McCLAIN 1/29/2020

PART NUMBER: AX069-A

DESCRIPTION: MAESTRO, DISSECTION BOARD, DARK BLUE W/ 3" LEGS

SIZE: A2	FILE NAME: AX069-A.dft	REV: A
SCALE: 1:3.03	DO NOT SCALE IF IN DOUBT ASK	SHEET 1 OF 2

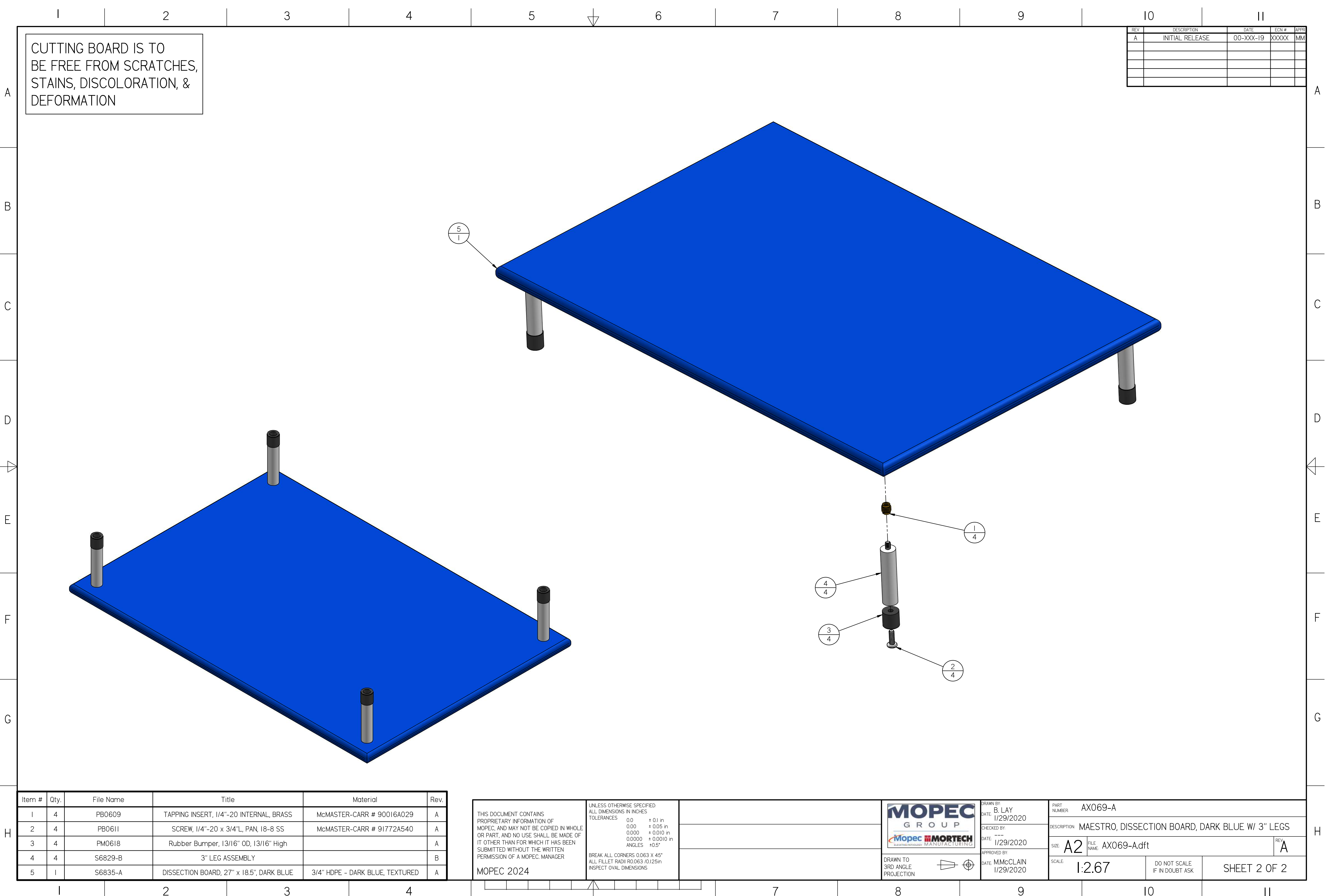

REV	DESCRIPTION	DATE	ECN #	APPR
A	INITIAL RELEASE	00-XXX-19	XXXXX	MM

Item #	Qty.	File Name	Title	Material	Rev.
1	4	PB0609	TAPPING INSERT, 1/4"-20 INTERNAL, BRASS	McMASTER-CARR # 90016A029	A
2	4	PB0611	SCREW, 1/4"-20 x 3/4"L, PAN, 18-8 SS	McMASTER-CARR # 91772A540	A
3	4	PM0618	Rubber Bumper, 13/16" OD, 13/16" High		A
4	4	S6829-B	3" LEG ASSEMBLY		B
5	1	S6835-A	DISSECTION BOARD, 27" x 18.5", DARK BLUE	3/4" HDPE - DARK BLUE, TEXTURED	A

THIS DOCUMENT CONTAINS  
PROPRIETARY INFORMATION OF  
MOPEC, AND MAY NOT BE COPIED IN WHOLE  
OR PART, AND NO USE SHALL BE MADE OF  
IT OTHER THAN FOR WHICH IT HAS BEEN  
SUBMITTED WITHOUT THE WRITTEN  
PERMISSION OF A MOPEC MANAGER

MOPEC 2024

UNLESS OTHERWISE SPECIFIED:  
ALL DIMENSIONS IN INCHES


TOLERANCES

0.0	± 0.1 in
0.00	± 0.05 in
0.000	± 0.010 in
0.0000	± 0.0010 in
ANGLES	± 0.5°

BREAK ALL CORNERS 0.063 X 45°  
ALL FILLET RADII R0.063 /0.125in  
INSPECT OVAL DIMENSIONS



DRAWN TO  
3RD ANGLE  
PROJECTION

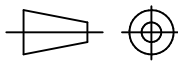

DRAWN BY:  
B. LAY  
DATE: 1/29/2020

CHECKED BY:  
DATE: 1/29/2020

APPROVED BY:  
DATE: M.McCLAIN  
1/29/2020

PART NUMBER:	AX069-A		
DESCRIPTION:	MAESTRO, DISSECTION BOARD, DARK BLUE W/ 3" LEGS		
SIZE:	A2	FILE NAME:	AX069-A.dft
SCALE:	1:2.67	DO NOT SCALE IF IN DOUBT ASK	SHEET 2 OF 2

\*AX069-A.dft\*