

REV	DESCRIPTION	DATE	ECN #	APPR
A	INITIAL RELEASE	00-XXX-19	XXXXX	MM

MOPEC LEGAL NOTE
THIS DOCUMENT CONTAINS
PROPRIETARY INFORMATION OF
MOPEC, AND MAY NOT BE COPIED IN WHOLE OR
PART, AND NO USE SHALL BE MADE OF IT
OTHER THAN FOR WHICH IT HAS BEEN
SUBMITTED WITHOUT THE WRITTEN
PERMISSION OF A MOPEC MANAGER
MOPEC 2024

UNLESS OTHERWISE SPECIFIED:
ALL DIMENSIONS IN INCHES
TOLERANCES
0.0 ± 0.1 in
0.00 ± 0.05 in
0.000 ± 0.010 in
0.0000 ± 0.0010 in
ANGLES ±0.5°
BREAK ALL CORNERS 0.063 X 45°
ALL FILLET RADII R0.063 /0.125in
INSPECT OVAL DIMENSIONS

MOPEC
GROUP

Mopec **MORTECH**
SCIENTIFIC TECHNOLOGY MANUFACTURING

DRAWN TO
3RD ANGLE
PROJECTION

DRAWN BY: jjohnson
DATE: 12/2/2024
CHECKED BY:
DATE: 12/2/2024
APPROVED BY:
DATE: M.McCLAIN
12/2/2024

PART NUMBER: HBI550-A	REV: A
DESCRIPTION: ELEVATING DISSECTION CART W/ LID ASSIST, ADA	
SIZE: A2	FILE NAME: HBI550-A.dft
SCALE: 1:8	DO NOT SCALE IF IN DOUBT ASK
	SHEET 1 OF 2



HB1550-A.dft

A
B
C
D
E
F
G
H

A
B
C
D
E
F
G
H

HOODS SHOWN IN CLOSED POSITION

HOODS SHOWN IN OPEN POSITION

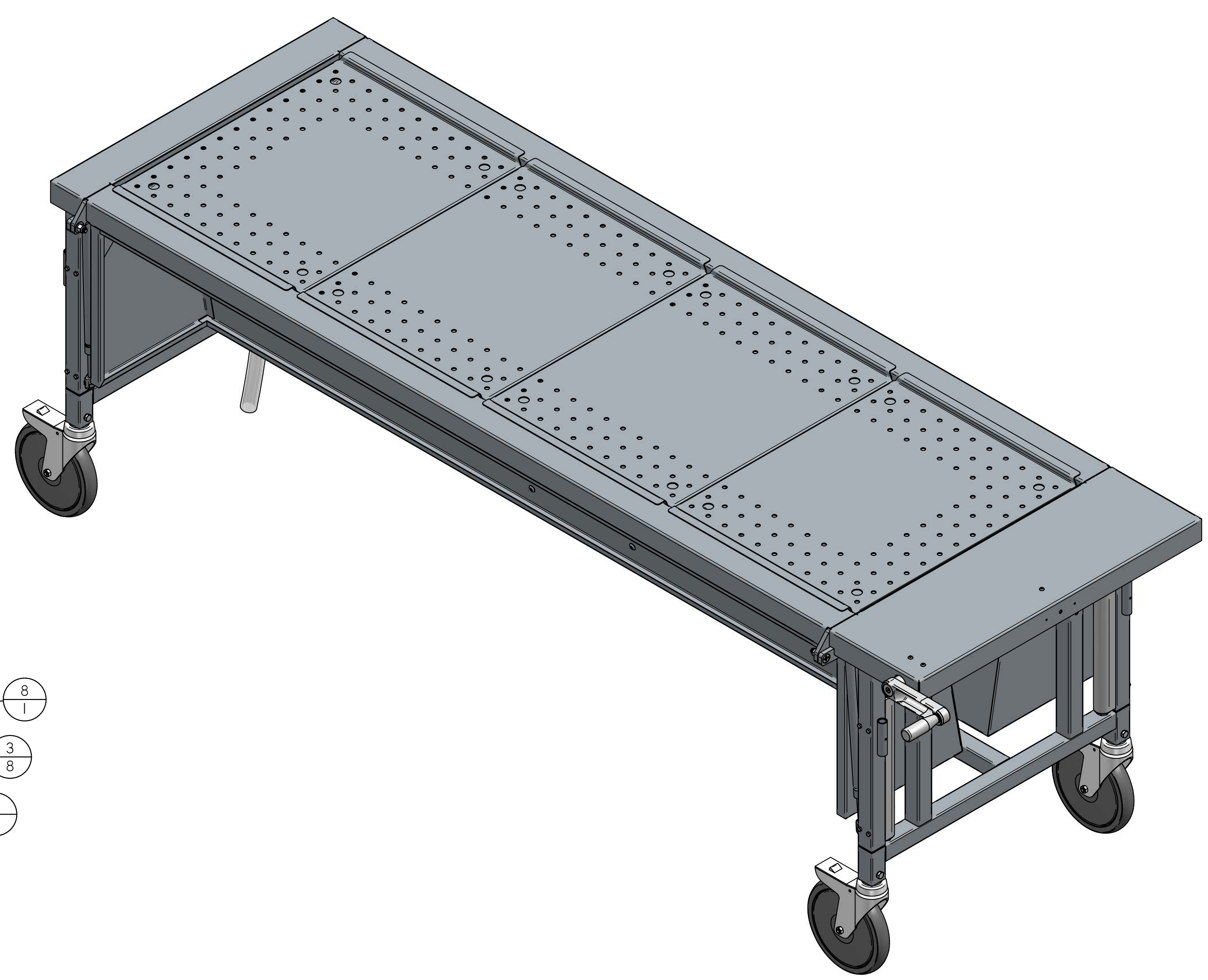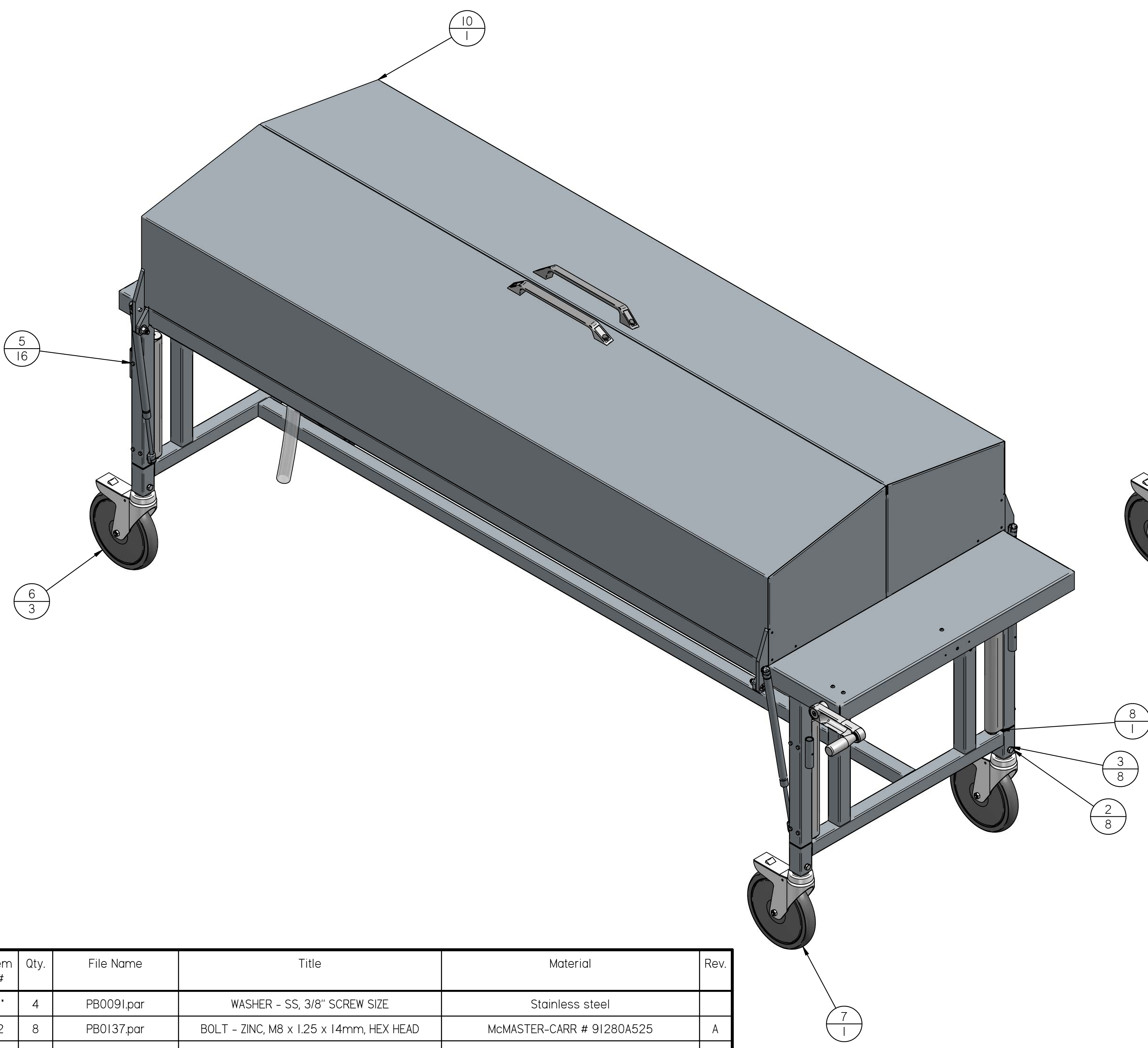

Item #	Qty.	File Name	Title	Material	Rev.
1*	4	PB009I.par	WASHER - SS, 3/8" SCREW SIZE	Stainless steel	
2	8	PB0137.par	BOLT - ZINC, M8 x 1.25 x 14mm, HEX HEAD	McMASTER-CARR # 91280A525	A
3	8	PB021I.par	WASHER - SS, M8 x 5/16", SERRATED BELLEVILLE	McMASTER-CARR # 90895A030	A
4*	4	PB0427.par	SCREW-CAP M10 X 65 MM FULLY THREADED SS	McMASTER-CARR # 91287A372	A
5	16	PB0603.par	BOLT HEX M5 X 45MM SS FULL THREAD	McMASTER-CARR # 91287A331	A
6	3	PD0026.par	CASTER - 8" TL (POWDER COATED) RED	TENTE # 2447UAP200R05-28	A
7	1	PD0027.par	CASTER - 8" DL (POWDER COATED) GREEN	TENTE # 2441UAP200R05-28	A
8	1	PY0006-A.asm	SUSPA, 200MM HC AND LIFTS		A
9*	4	S2290-E.par	CASTER ADAPTER, 1-1/2" SQ. 16 GA. TUBE	RM0620, TUBING, RND, MS, 1-3/8" OD X 10 GA, DOM	E
10	1	S6940-A.asmICLOSED	HBI500 ELEVATING DISSECTION CART ASSEMBLY		A

MOPEC LEGAL NOTE
THIS DOCUMENT CONTAINS PROPRIETARY INFORMATION OF MOPEC, AND MAY NOT BE COPIED IN WHOLE OR PART, AND NO USE SHALL BE MADE OF IT OTHER THAN FOR WHICH IT HAS BEEN SUBMITTED WITHOUT THE WRITTEN PERMISSION OF A MOPEC MANAGER
MOPEC 2024

UNLESS OTHERWISE SPECIFIED:
ALL DIMENSIONS IN INCHES
TOLERANCES
0.0 ± 0.1 in
0.00 ± 0.05 in
0.000 ± 0.010 in
0.0000 ± 0.0010 in
ANGLES ± 0.5°
BREAK ALL CORNERS 0.063 X 45°
ALL FILLET RADII R0.063 (0.125 in)
INSPECT OVAL DIMENSIONS



DRAWN BY: jjohnson
DATE: 12/2/2024
CHECKED BY:
DATE: 12/2/2024
APPROVED BY:
DATE: M.McCLAIN 12/2/2024

PART NUMBER: HBI550-A	DESCRIPTION: ELEVATING DISSECTION CART W/ LID ASSIST, ADA	REV: A
SIZE: A2	FILE NAME: HBI550-A.dft	
SCALE: 1:8	DO NOT SCALE IF IN DOUBT ASK	SHEET 2 OF 2

HB1550-A.dft