

REV	DESCRIPTION	DATE	ECN #	APPR
A	INITIAL RELEASE	00-XXX-19	XXXXX	MM

Item Number	File Name (no extension)	Author	Quantity	Material
1	PB0609	jbourbeau	4	McMASTER-CARR # 90016A029
2	S6835-A	jbourbeau	1	3/4" HDPE - DARK BLUE, TEXTURED

THIS DOCUMENT CONTAINS PROPRIETARY INFORMATION OF MOPEC, AND MAY NOT BE COPIED IN WHOLE OR PART, AND NO USE SHALL BE MADE OF IT OTHER THAN FOR WHICH IT HAS BEEN SUBMITTED WITHOUT THE WRITTEN PERMISSION OF A MOPEC MANAGER


MOPEC 2024

UNLESS OTHERWISE SPECIFIED:
ALL DIMENSIONS IN INCHES

TOLERANCES

0.0	± 0.1 in
0.00	± 0.05 in
0.000	± 0.010 in
0.0000	± 0.0010 in
ANGLES	± 0.5°


BREAK ALL CORNERS 0.063 X 45°
ALL FILLET RADII R0.063 (0.125in)
INSPECT OVAL DIMENSIONS



MOPEC GROUP

subcontractor MORTTECH MANUFACTURING

DRAWN TO 3RD ANGLE PROJECTION

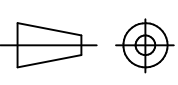

DRAWN BY: B. LAY
DATE: 1/30/2025

CHECKED BY: _____
DATE: 1/30/2025

APPROVED BY: _____
DATE: M.McCLAIN 1/30/2025

PART NUMBER: AX068-060-A

DESCRIPTION: MAESTRO, DISSECTION BOARD, DARK BLUE

SIZE: A2	FILE NAME: AX068-060-A.dft	REV: A
SCALE: 1:2.67	DO NOT SCALE IF IN DOUBT ASK	SHEET 2 OF 2