
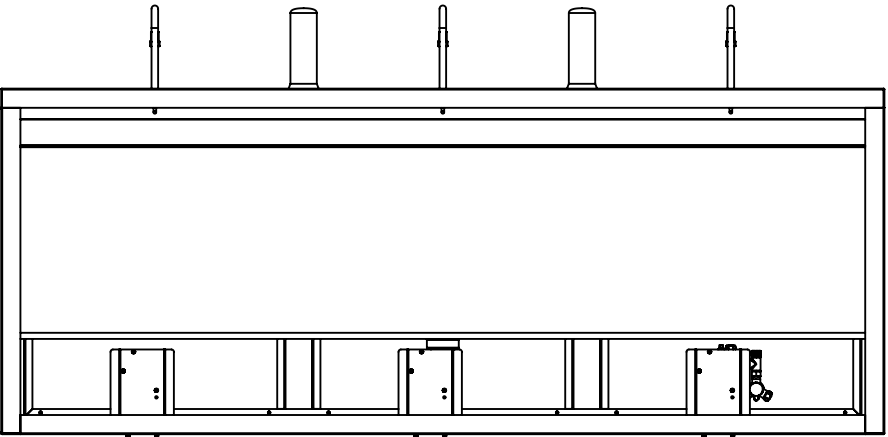
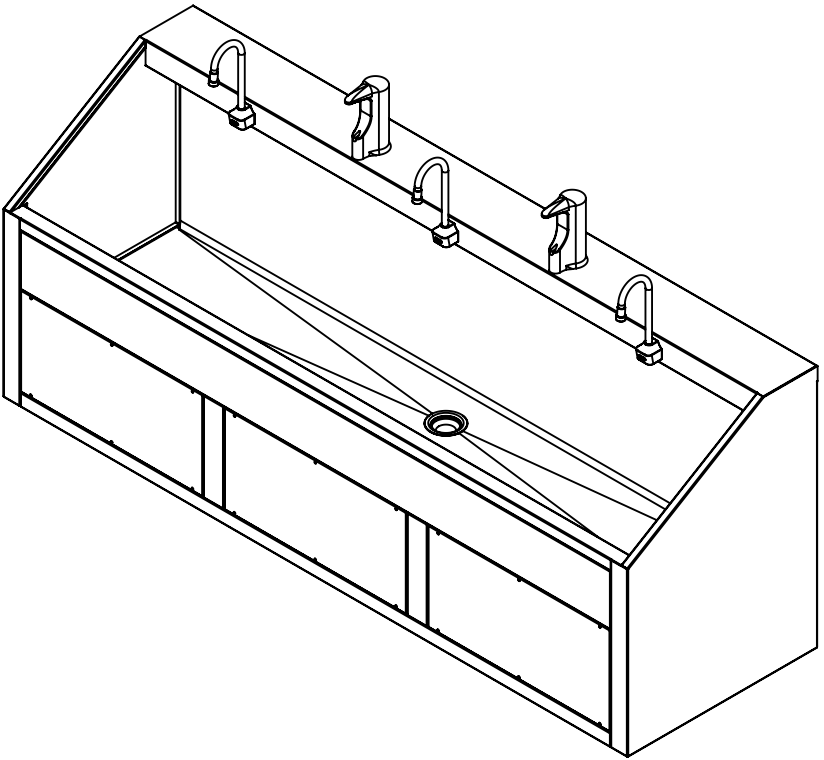
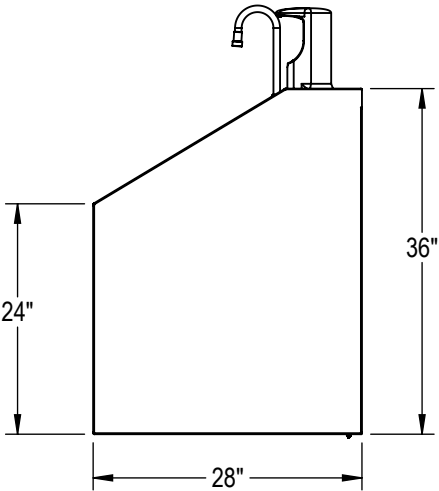
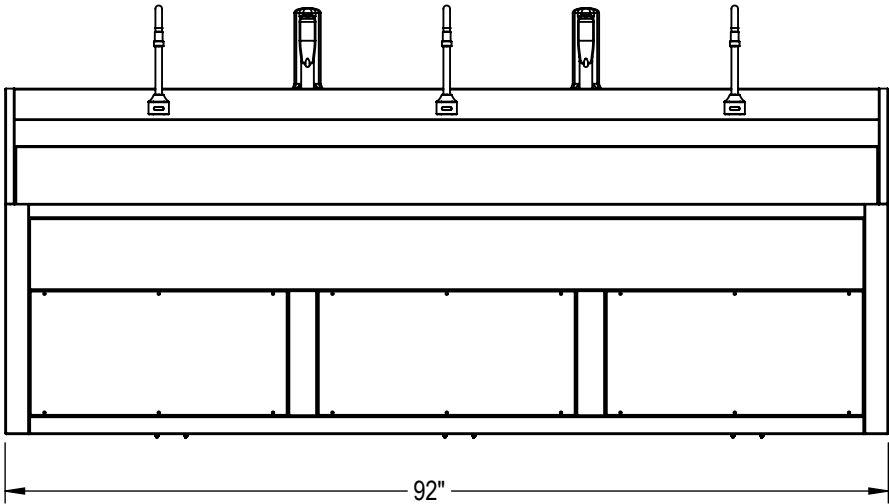


NOTES: UNLESS OTHERWISE SPECIFIED


1. APPLICABLE TO STANDARDS AND SPECIFICATIONS;
ASME Y14.5-2009, DIMENSIONING AND TOLERANCING
ASME Y14.38-2007, ABBREVIATIONS

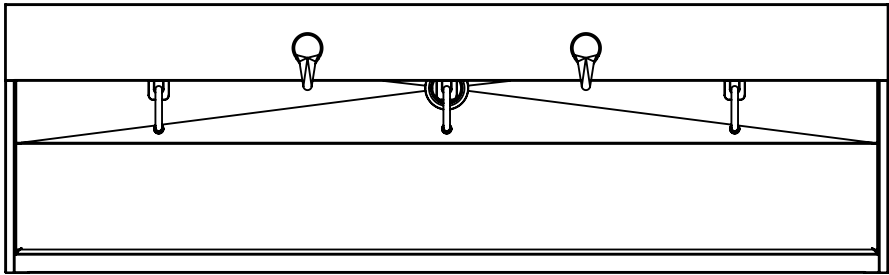

TOLERANCE BLOCK:		PROPRIETARY THIS DOCUMENT AND THE DATE DISCLOSED HEREIN AND HEREWITH IS NOT TO BE REPRODUCED, USED, OR OTHERWISE DISCLOSED IN WHOLE OR IN PART TO ANYONE WITHOUT WRITTEN AUTHORIZATION FROM MORTECH MFG. MOPEC GROUP Mopec ELEVATING PATHOLOGY MORTECH MANUFACTURING		TITLE							
DIMENSIONS ARE IN INCHES											
DECIMALS .X ± .1	ANGLES ± 0° 30'										
DECIMALS .XX ± .06	FRACTIONS ± 1/16"										
DECIMALS .XXX ± .031		APPROVALS		DATE		SIZE B DWG NO. SS103 REV. X01					
DO NOT SCALE DRAWING		DRAWN BY									
MATERIAL		CHECKED BY									
304 STAINLESS STEEL											
FINISH		APPROVED BY				SCALE		NEXT ASSY.		SHEET	
#4 SATIN						1:20				1 OF 2	

NOTES: UNLESS OTHERWISE SPECIFIED

1. APPLICABLE TO STANDARDS AND SPECIFICATIONS;
ASME Y14.5-2009, DIMENSIONING AND TOLERANCING
ASME Y14.38-2007, ABBREVIATIONS

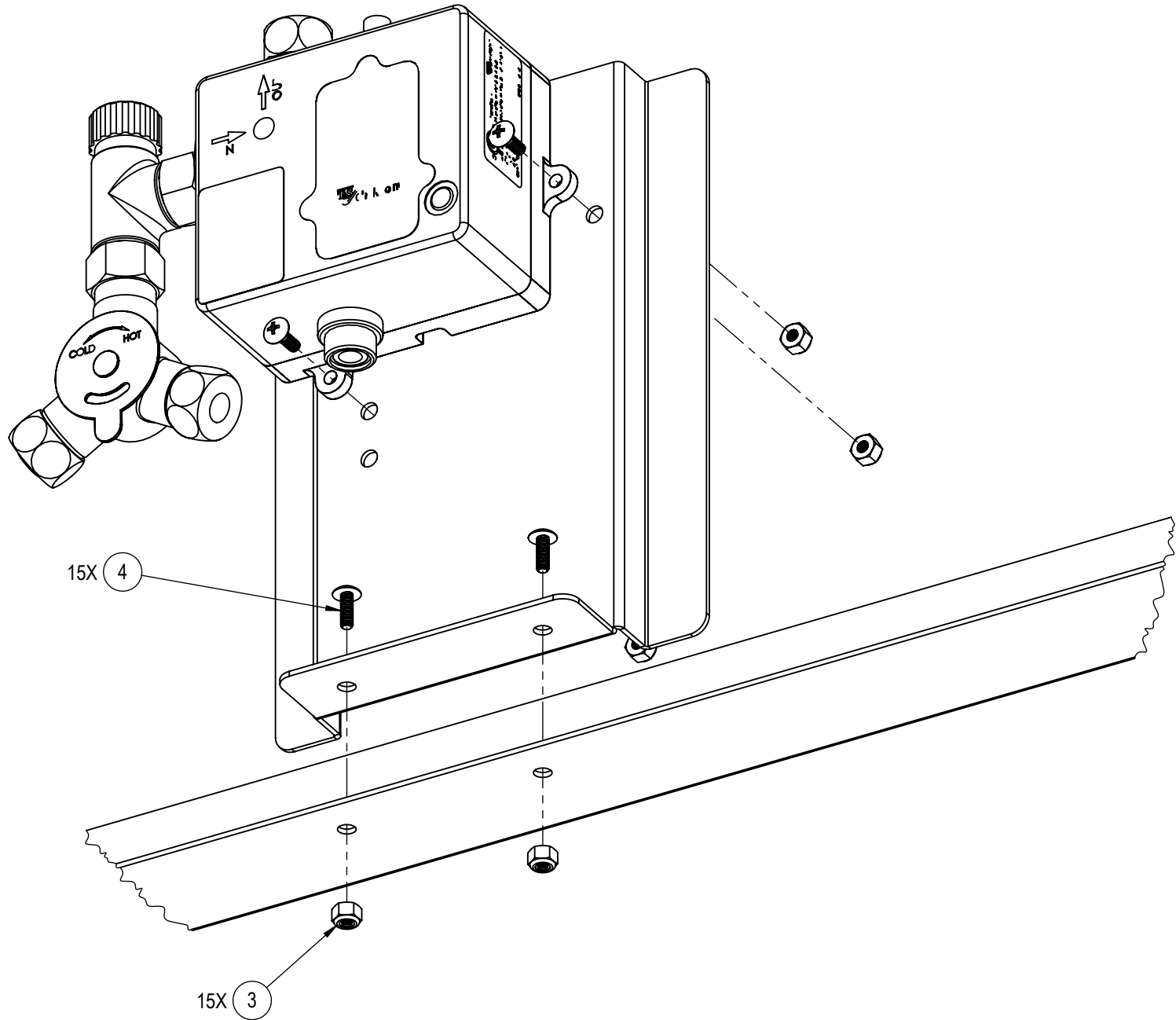

9	3	SS101-ACCESS-PANEL	PANEL, ACCESS- 13.00 X 26.81
8	1	STRAINERS350	STRAINER, SINK- FOR 3.50" OPENING
7	2	DISPENSERSOAP	
6	3	FAUCETWM11	FAUCET, WALL MOUNT- 11" X 4 1/8"
5	1	EC3101_CONTROL	
4	15	SCREWPTH8325SS	SCREW, PHILLIPS TRUSS HEAD- #8-32 X .500 L, SS
3	15	NUTLK832S	NUT, LOCK- 8-32, WITH NYLON INSERT, SS
2	3	SSHF-8000	
1	1	SS103-7000	TRIPLE SCRUB STATION, W/ HANDS FREE, SUBASSEMBLY
ITEM NO.	QTY.	PART NUMBER	DESCRIPTION

PARTS LIST

TOLERANCE BLOCK: DIMENSIONS ARE IN INCHES		PROPRIETARY THIS DOCUMENT AND THE DATE DISCLOSED HEREIN AND HERewith IS NOT TO BE REPRODUCED, USED, OR OTHERWISE DISCLOSED IN WHOLE OR IN PART TO ANYONE WITHOUT WRITTEN AUTHORIZATION FROM MORTECH MFG.		MOPEC GROUPMopecMORTECH ELEVATING PATHOLOGY MANUFACTURING	
DECIMALS .X ± .1	ANGLES ± 0° 30'	APPROVALS		DATE	
DECIMALS .XX ± .06	FRACTIONS ± 1/16"	DO NOT SCALE DRAWING		DRAWN BY	
DECIMALS .XXX ± .031		304 STAINLESS STEEL		CHECKED BY	
FINISH	#4 SATIN	APPROVED BY		SIZE B	
				DWG NO. SS103	
				REV. X01	
				SCALE 1:2	
				NEXT ASSY.	
				SHEET 2 OF 2	