


PARTS ARE FOR KITTING PURPOSE ONLY

REV	DESCRIPTION	DATE	ECN #	APPR
A	INITIAL RELEASE	00-XXX-19	XXXXX	MM

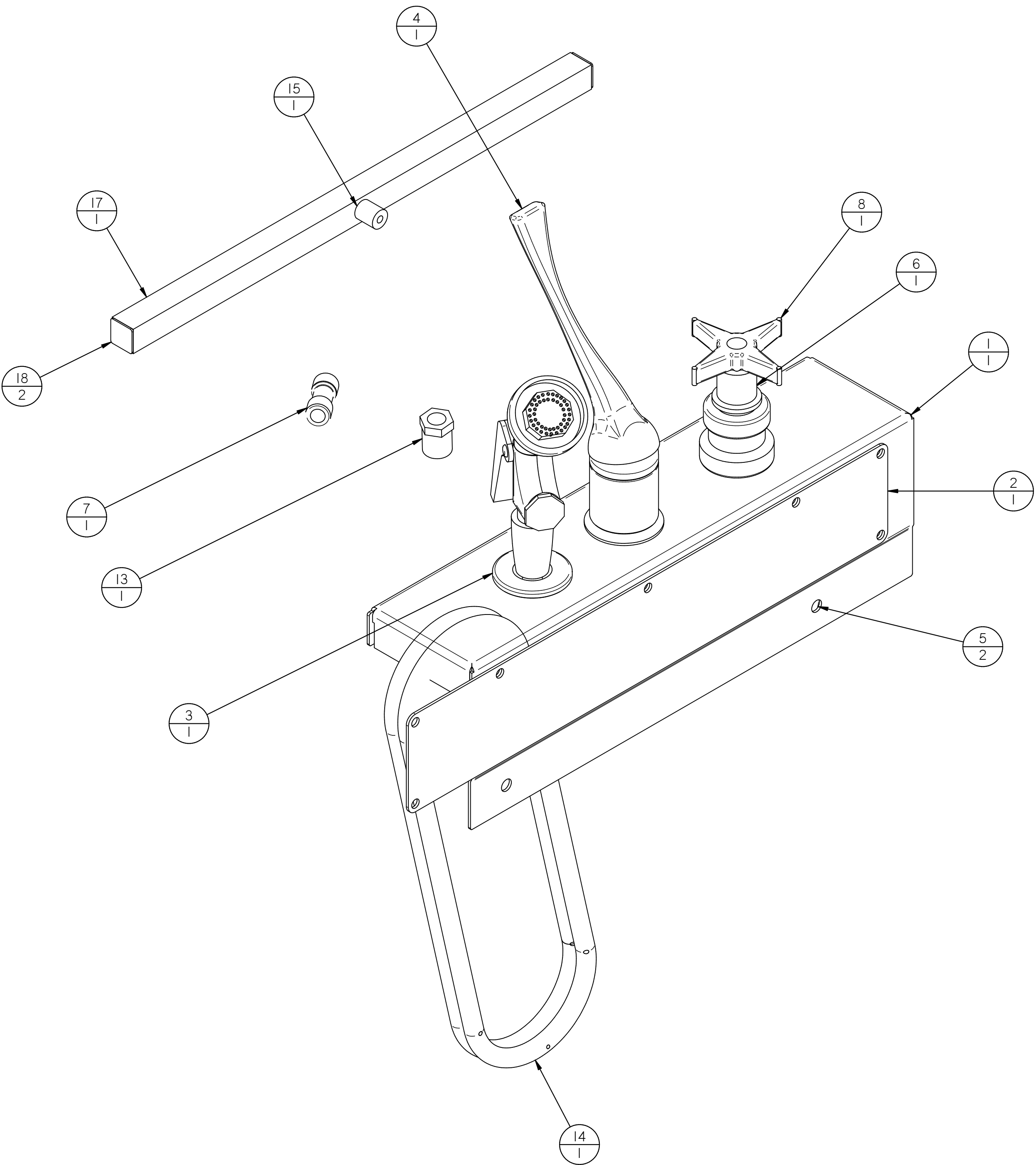

Item #	Qty.	File Name	Title	Material	Rev.
1	1	M-5111-B.psm	MB, WORK SURFACE, RIGHT END CAP	14 GA. TYPE T304 SS	B
2	1	M-5113-B.psm	MB,END CAP COVER	SHT, SS, 14 GA, 304	B
3	1	PP0056.par	FIXTURE - BRASS DECK FLANGE - 1" ID. X 3" L	FISHER #F00213	A
4	1	PP0079.par	FIXTURE DECK MTD. H/C SGL. LEVER		A
5	2	PP0199.par	FITTING BRASS B/F PREVENTER 3/8" FNPT		A
6	1	PP0654.par	WATER SAVER VALVE	WATER SAVER	A
7	1	PP0694.par	UNION TEE, TUBE 3/8	PARKER #6304 60 00WP2	A
8	1	PP0941.par	STAR HANDLE FOR VALVE	WATER SAVER	A
9*	3	PPI011.par	UNION TEE, TUBE 3/8	PARKER #6304 60 00WP2	A
10*	2	PPI013.par	UNION ELBOW, TUBE 3/8	PARKER #6302 60 00WP2	A
11*	2	PPI014.par	MALE, NPTF THD, TUBE 3/8- 3/8 NPTF	PARKER #6505 60 18WP2	A
12*	2	PPI016.par	SWIVEL ELBOW NPTF, TUBE 3/8-3/8	PARKER #6509 60 18 WP2	A
13	1	PPI035.par	Bushing Reducing Adapter, 1/2 Male x 1/4 Female NPTF	MCMaster CARR : #9151K95	A
14	1	PPI038.asm	HOSEASSY1.asm		A
15	1	PPI046.par	Fully Threaded, 1/4 NPT, 7/8" Long	MCMaster CARR: 4830K131	A
16*	1	PPI078.par			
17	1	S6848-A.par	MB, END RINSE BAR	RM0624, TUBING , SQ, SS, 3/4" X 16 GA, 304 , P	A
18	2	S6849-A.psm	MB, END RINSE CAP	18 GA. TYPE T304 SS	A

THIS DOCUMENT CONTAINS PROPRIETARY INFORMATION OF MOPEC, AND MAY NOT BE COPIED IN WHOLE OR PART, AND NO USE SHALL BE MADE OF IT OTHER THAN FOR WHICH IT HAS BEEN SUBMITTED WITHOUT THE WRITTEN PERMISSION OF A MOPEC MANAGER MOPEC 2020	UNLESS OTHERWISE SPECIFIED: ALL DIMENSIONS IN INCHES TOLERANCES 0.0 ± 0.1 in 0.00 ± 0.05 in 0.000 ± 0.010 in 0.0000 ± 0.0010 in ANGLES ±0.5° BREAK ALL CORNERS 0.063 X 45° ALL FILLET RADII R0.063 /0.125in INSPECT OVAL DIMENSIONS	MATERIAL: MATERIAL USAGE: Error: No reference	 	DRAWN BY: DATE: jjohnson 4/15/2025	PART NUMBER: MB1036L-A		
		FINISH: Error: No reference		CHECKED BY: DATE: --- 4/15/2025	DESCRIPTION: MAESTRO, HANDSPRAY & END RINSE, LEFT HAND	SIZE: A2 FILE NAME: MB1036L-A.dft	REV: A
				DRAWN TO 3RD ANGLE PROJECTION 	APPROVED BY: DATE: MMcCLAIN 4/15/2025	SCALE: 1:2	DO NOT SCALE IF IN DOUBT ASK

MB1036L-A.dft