


PARTS ARE FOR KITTING PURPOSE ONLY

REV	DESCRIPTION	DATE	ECN #	APPR
A	INITIAL RELEASE	00-XXX-19	XXXXX	MM

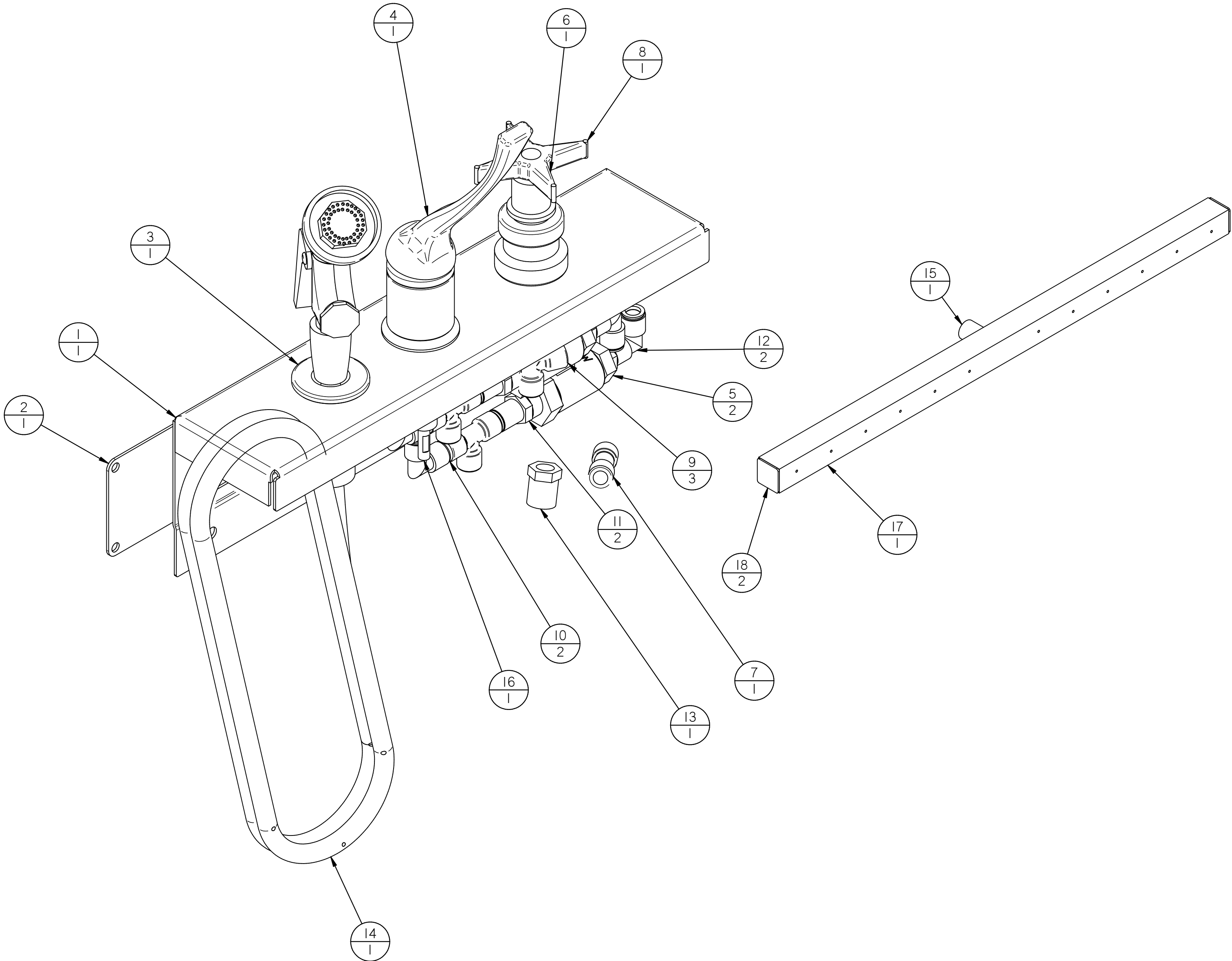

Item #	Qty.	File Name	Title	Material	Rev.
1	1	M-5111-02-B	MB, WORK SURFACE, LEFT END CAP	SHT, SS, 14 GA, 304	B
2	1	M-5113-B	MBEND CAP COVER	SHT, SS, 14 GA, 304	B
3	1	PP0056	FITURE - BRASS DECK FLANGE - 1" ID. X 3" L	FISHER #F00213	A
4	1	PP0079	FITURE DECK MTD. H/C SGL. LEVER		A
5	2	PP0199	FITTING BRASS B/F PREVENTER 3/8" FNPT		A
6	1	PP0654	WATER SAVER VALVE	WATER SAVER	A
7	1	PP0694	UNION TEE, TUBE 3/8	PARKER #6304 60 00WP2	A
8	1	PP0941	STAR HANDLE FOR VALVE	WATER SAVER	A
9	3	PPI011	UNION TEE, TUBE 3/8	PARKER #6304 60 00WP2	A
10	2	PPI013	UNION ELBOW, TUBE 3/8	PARKER #6302 60 00WP2	A
11	2	PPI014	MALE, NPTF THD, TUBE 3/8- 3/8 NPTF	PARKER #6505 60 18WP2	A
12	2	PPI016	SWIVEL ELBOW NPTF, TUBE 3/8-3/8	PARKER #6509 60 18 WP2	A
13	1	PPI035	Bushing Reducing Adapter, 1/2 Male x 1/4 Female NPTF	MCMaster CARR : #9151K95	A
14	1	PPI038	HOSEASSY1.asm		A
15	1	PPI046	Fully Threaded, 1/4 NPT, 7/8" Long	MCMaster CARR: 4830K131	A
16	1	PPI078			
17	1	S6848-A	MB, END RINSE BAR	RM0624, TUBING , SQ, SS, 3/4" X 16 GA, 304 , P	A
18	2	S6849-A	MB, END RINSE CAP	18 GA. TYPE T304 SS	A

THIS DOCUMENT CONTAINS PROPRIETARY INFORMATION OF MOPEC, AND MAY NOT BE COPIED IN WHOLE OR PART, AND NO USE SHALL BE MADE OF IT OTHER THAN FOR WHICH IT HAS BEEN SUBMITTED WITHOUT THE WRITTEN PERMISSION OF A MOPEC. MANAGER

MOPEC 2020

UNLESS OTHERWISE SPECIFIED:
ALL DIMENSIONS IN INCHES
TOLERANCES
0.0 ± 0.1 in
0.00 ± 0.05 in
0.000 ± 0.010 in
0.0000 ± 0.0010 in
ANGLES ±0.5°
BREAK ALL CORNERS 0.063 X 45°
ALL FILLET RADII R0.063 /0.125in
INSPECT OVAL DIMENSIONS

MATERIAL:
MATERIAL USAGE:
Error: No reference
FINISH:
Error: No reference

MOPEC GROUP

MOPEC MORTTECH

DRAWN TO 3RD ANGLE PROJECTION

DRAWN BY: jjohnson

DATE: 4/15/2025

CHECKED BY:

DATE: 4/15/2025

APPROVED BY:

DATE: M.McCLAIN 4/15/2025

PART NUMBER: MB1036R-A

DESCRIPTION: MAESTRO, HANDSPRAY & END RINSE, RIGHT HAND

SIZE: A2

FILE NAME: MB1036R-Adft

SCALE: 1:2

DO NOT SCALE IF IN DOUBT ASK

SHEET 1 OF 1

MB1036R-A.dft