



MB1039-096-B.dft

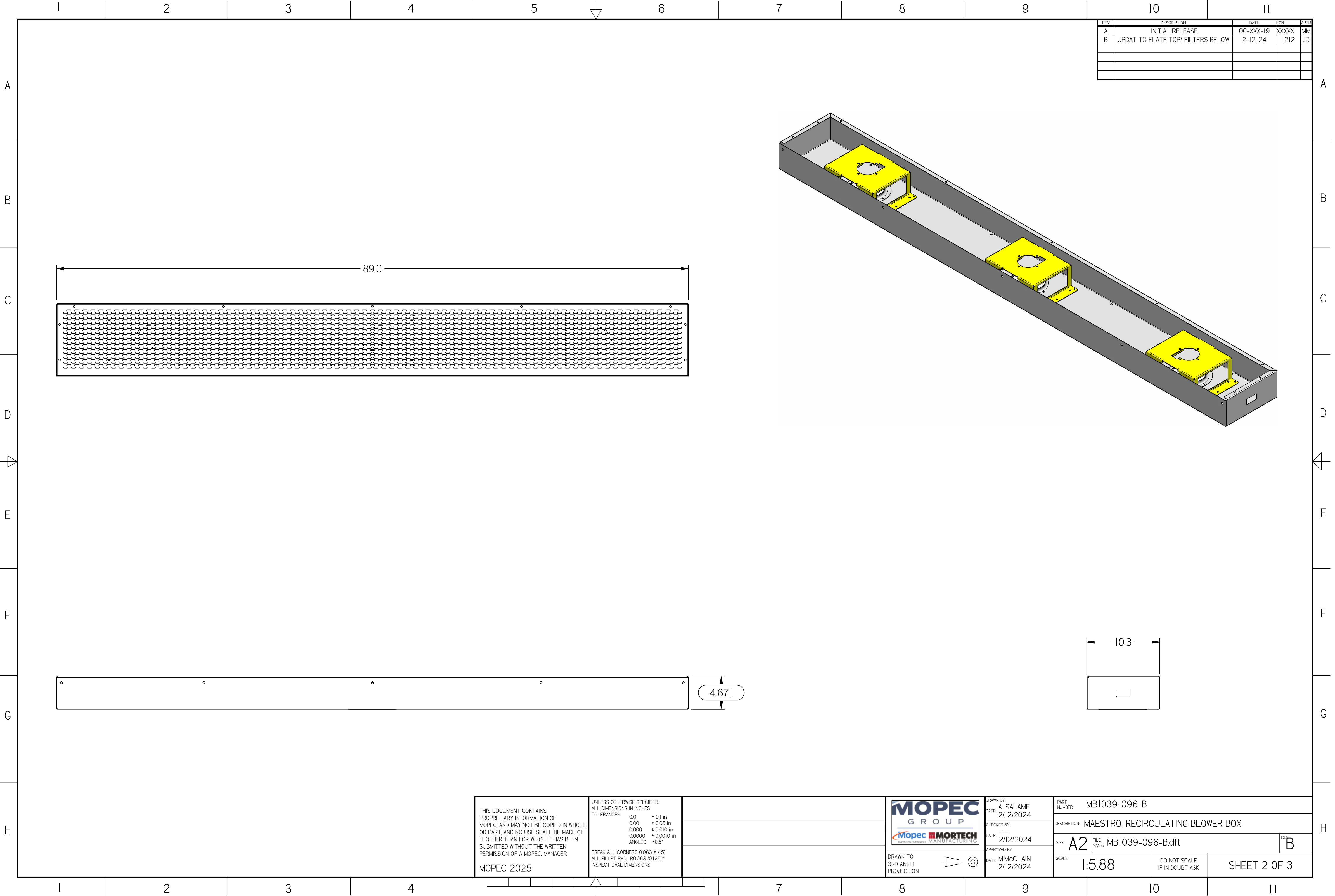

MB1039-096-B.dft

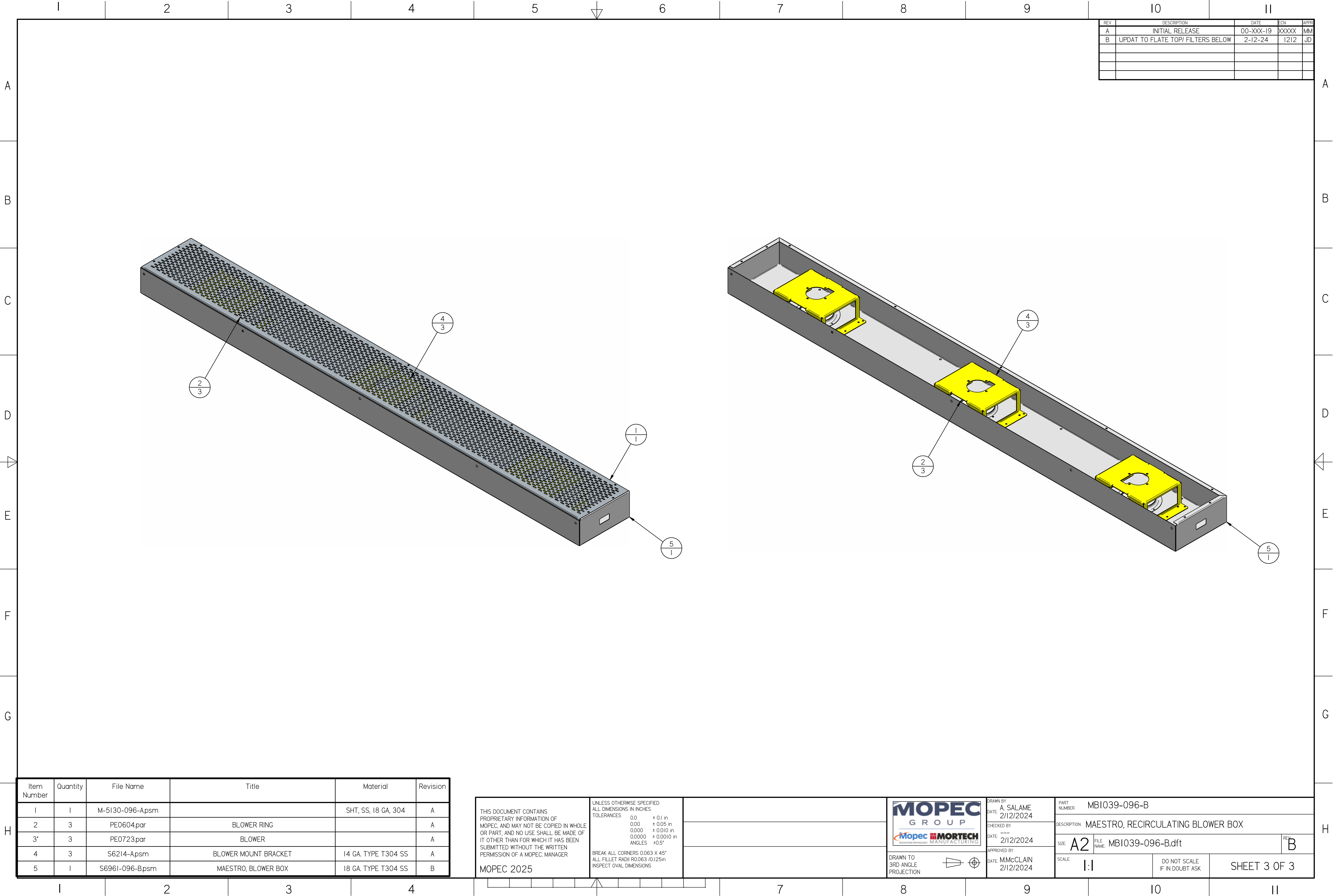

REV	DESCRIPTION	DATE	ECN	APPR
A	INITIAL RELEASE	00-XXX-19	XXXXX	MM
B	UPDAT TO FLATE TOP/ FILTERS BELOW	2-12-24	I212	JD

Item Number	Quantity	File Name	Title	Material	Revision
1	1	M-5130-096-Apsm		SHT, SS, 18 GA, 304	A
2	3	PE0604.par	BLOWER RING		A
3	3	PE0723.par	BLOWER		A
4	3	S6214-Apsm	BLOWER MOUNT BRACKET	14 GA. TYPE T304 SS	A
5	1	S6961-096-B.psm	MAESTRO, BLOWER BOX	18 GA. TYPE T304 SS	B

THIS DOCUMENT CONTAINS
PROPRIETARY INFORMATION OF
MOPEC, AND MAY NOT BE COPIED IN WHOLE
OR PART, AND NO USE SHALL BE MADE OF
IT OTHER THAN FOR WHICH IT HAS BEEN
SUBMITTED WITHOUT THE WRITTEN
PERMISSION OF A MOPEC. MANAGER

MOPEC 2025

UNLESS OTHERWISE SPECIFIED:
ALL DIMENSIONS IN INCHES


TOLERANCES

0.0	± 0.1 in
0.00	± 0.05 in
0.000	± 0.010 in
0.0000	± 0.0010 in
ANGLES	± 0.5°

BREAK ALL CORNERS 0.063 X 45°
ALL FILLET RADII R0.063 /0.125in
INSPECT OVAL DIMENSIONS



DRAWN TO
3RD ANGLE
PROJECTION

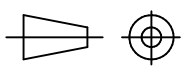

DRAWN BY:
A. SALAME
DATE: 2/12/2024

CHECKED BY:
DATE: ---
2/12/2024

APPROVED BY:
DATE: M.McCLAIN
2/12/2024

PART NUMBER:	MB1039-096-B		
DESCRIPTION:	MAESTRO, RECIRCULATING BLOWER BOX		
SIZE:	A2	FILE NAME:	MB1039-096-B.dft
SCALE:	1:1	DO NOT SCALE IF IN DOUBT ASK	SHEET 3 OF 3

MB1039-096-B.dft