



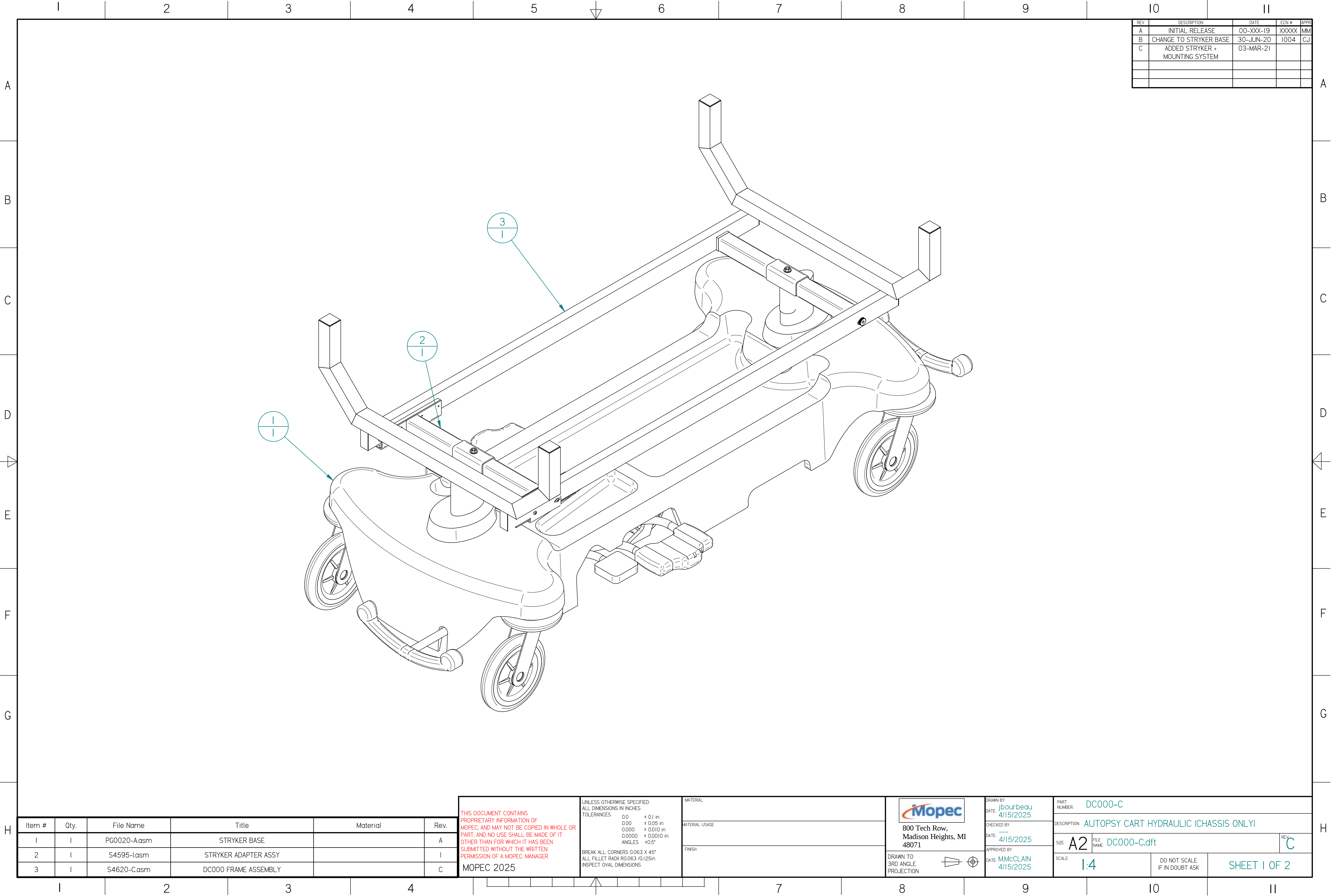

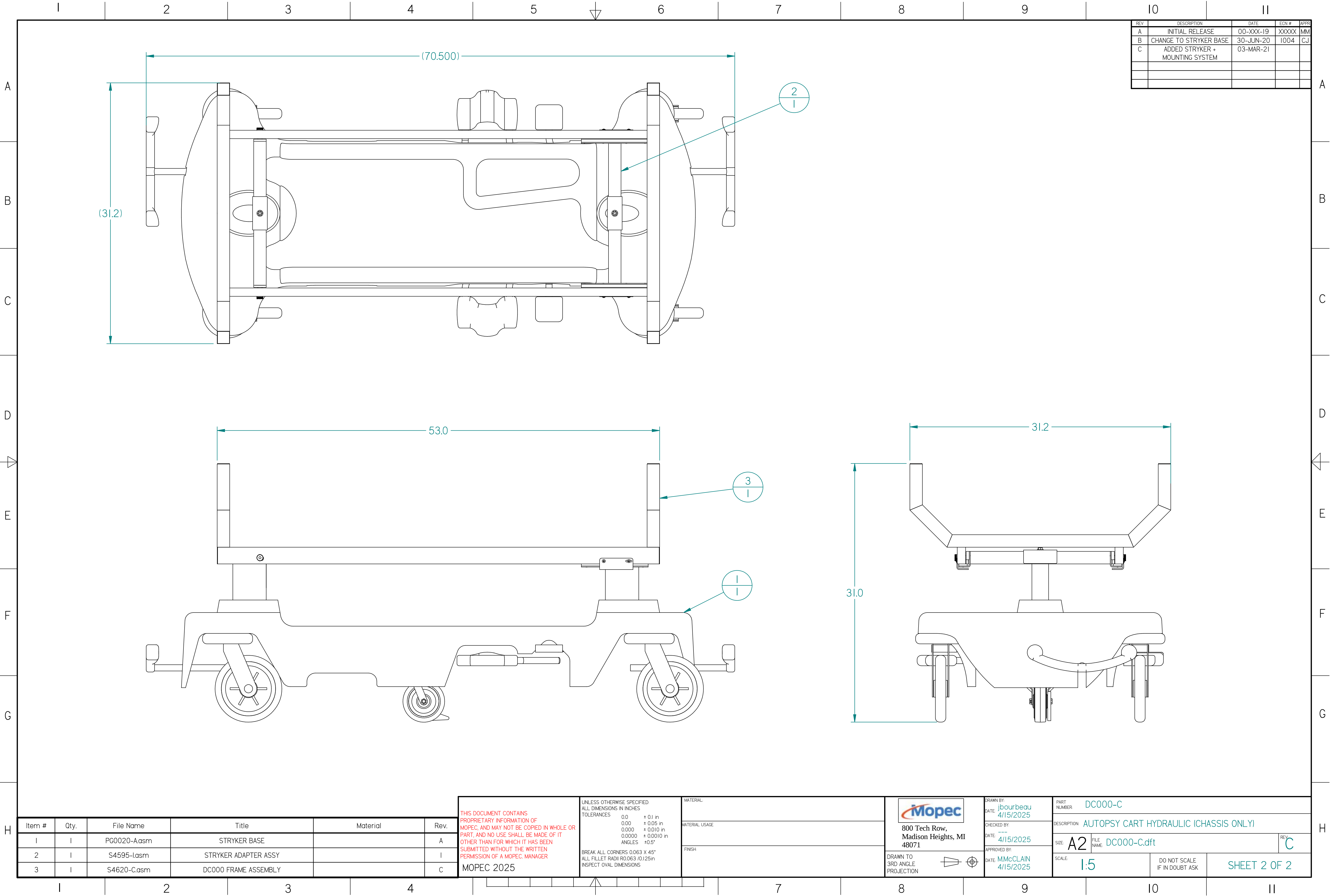

REV	DESCRIPTION	DATE	ECN #	APPR
A	INITIAL RELEASE	00-XXX-19	XXXXX	MM
B	CHANGE TO STRYKER BASE	30-JUN-20	1004	CJ
C	ADDED STRYKER + MOUNTING SYSTEM	03-MAR-21		

Item #	Qty.	File Name	Title	Material	Rev.
1	1	PG0020-A.asm	STRYKER BASE		A
2	1	S4595-I.asm	STRYKER ADAPTER ASSY		I
3	1	S4620-C.asm	DC000 FRAME ASSEMBLY		C

THIS DOCUMENT CONTAINS PROPRIETARY INFORMATION OF MOPEC, AND MAY NOT BE COPIED IN WHOLE OR PART, AND NO USE SHALL BE MADE OF IT OTHER THAN FOR WHICH IT HAS BEEN SUBMITTED WITHOUT THE WRITTEN PERMISSION OF A MOPEC MANAGER

MOPEC 2025

UNLESS OTHERWISE SPECIFIED:
ALL DIMENSIONS IN INCHES

TOLERANCES

0.0	± 0.1 in
0.00	± 0.05 in
0.000	± 0.010 in
0.0000	± 0.0010 in
ANGLES	± 0.5°

BREAK ALL CORNERS 0.063 X 45°
ALL FILLET RADII R0.063 / 0.125 in
INSPECT OVAL DIMENSIONS

MATERIAL

MATERIAL USAGE

FINISH

800 Tech Row,
Madison Heights, MI
48071

DRAWN TO 3RD ANGLE PROJECTION

DRAWN BY: jbourbeau
DATE: 4/15/2025

CHECKED BY: ---
DATE: 4/15/2025

APPROVED BY: MMcCLAIN
DATE: 4/15/2025

PART NUMBER: DC000-C

DESCRIPTION: AUTOPSY CART HYDRAULIC (CHASSIS ONLY)

SIZE: A2

SCALE: 1:5

DO NOT SCALE IF IN DOUBT ASK

SHEET 2 OF 2