
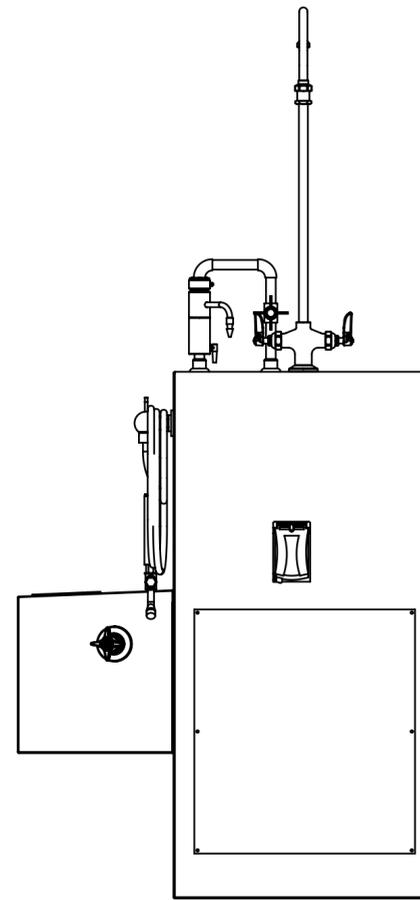
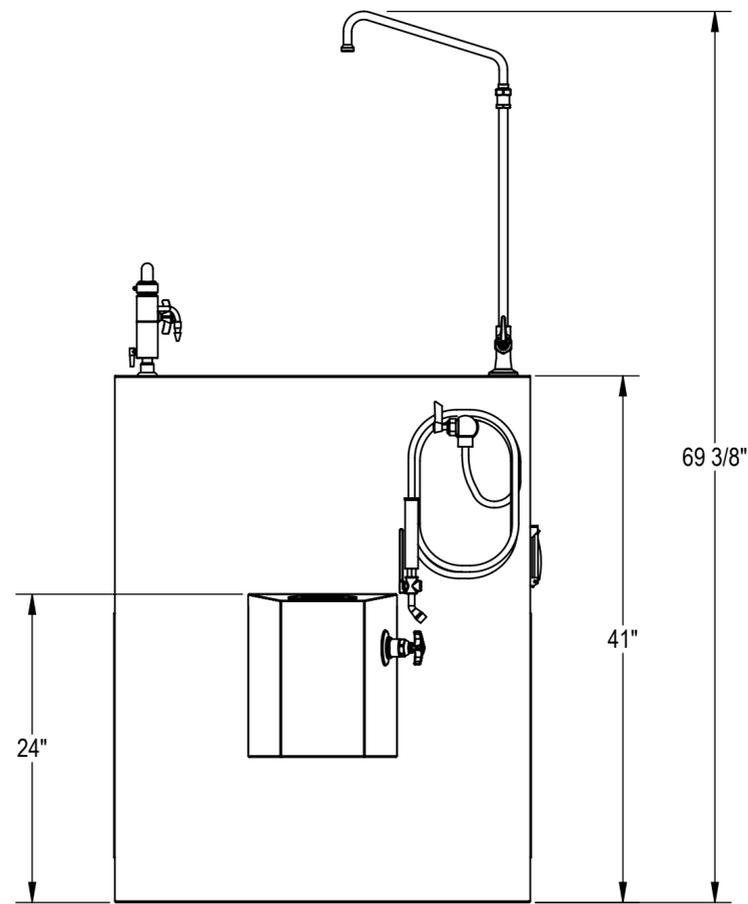
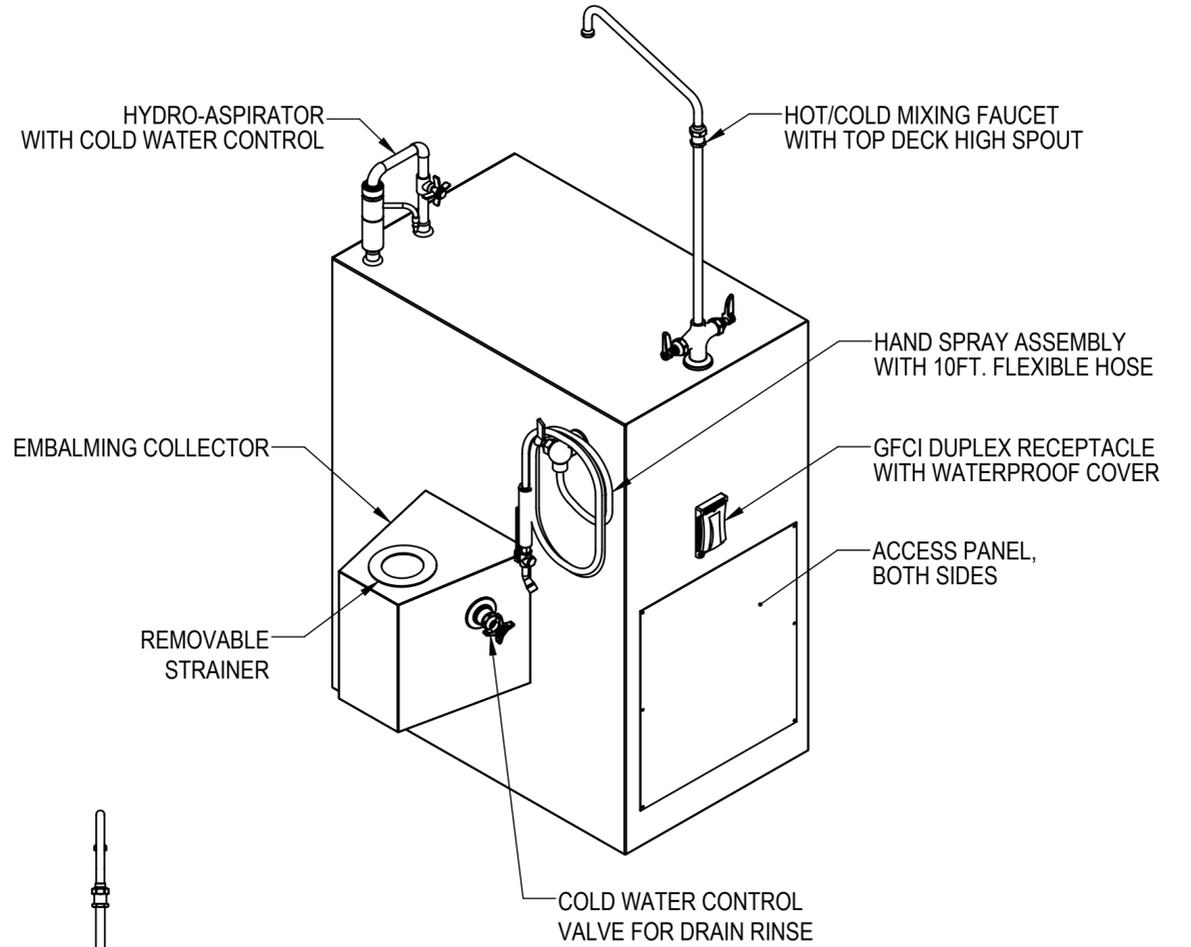


NOTES: UNLESS OTHERWISE SPECIFIED


1. APPLICABLE STANDARDS AND SPECIFICATIONS;
ANSI Y14.5M, DIMENSIONS AND TOLERANCES
ANSI Y1.1, ABBREVIATIONS

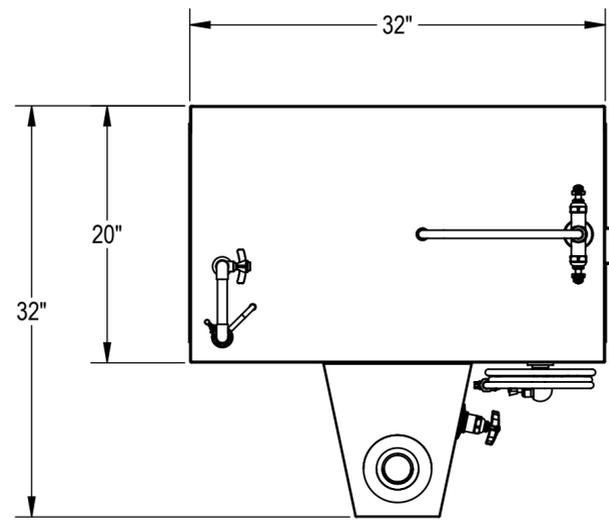

CUSTOMER		APPROVAL		<p style="text-align: center;">PROPRIETARY</p> <p style="text-align: center;">THIS DOCUMENT AND THE DATE DISCLOSED HEREIN AND HEREWITH IS NOT TO BE REPRODUCED, USED, OR OTHERWISE DISCLOSED IN WHOLE OR IN PART TO ANYONE WITHOUT WRITTEN AUTHORIZATION FROM MORTECH MFG.</p>			<p style="font-size: small;">411 North Aerojet Avenue Azusa, CA 91702 tel: (800) 410-0100 (626) 334-1471 fax: (626) 334-1704 www.mortechmfg.com</p>												
PROJECT									<p style="text-align: center;">DO NOT SCALE DRAWING</p> <table border="1" style="width: 100%; border-collapse: collapse;"> <tr> <th style="width: 33%;">MATERIAL</th> <th style="width: 33%;">APPROVALS</th> <th style="width: 33%;">DATE</th> </tr> <tr> <td>304 STAINLESS STEEL</td> <td>DRAWN BY GABRIEL G.</td> <td>03/10/17</td> </tr> <tr> <td>FINISH #4 SATIN</td> <td>CHECKED BY</td> <td></td> </tr> <tr> <td></td> <td>APPROVED BY</td> <td></td> </tr> </table>		MATERIAL	APPROVALS	DATE	304 STAINLESS STEEL	DRAWN BY GABRIEL G.	03/10/17	FINISH #4 SATIN	CHECKED BY	
MATERIAL	APPROVALS	DATE																	
304 STAINLESS STEEL	DRAWN BY GABRIEL G.	03/10/17																	
FINISH #4 SATIN	CHECKED BY																		
	APPROVED BY																		
		SIGNATURE		DATE		<table border="1" style="width: 100%; border-collapse: collapse;"> <tr> <th style="width: 25%;">SIZE</th> <th style="width: 25%;">DWG. NO.</th> <th style="width: 25%;">REV.</th> <th style="width: 25%;">REV.</th> </tr> <tr> <td>C</td> <td>10363-AP</td> <td></td> <td>X01</td> </tr> </table>		SIZE	DWG. NO.	REV.	REV.	C	10363-AP		X01				
SIZE	DWG. NO.	REV.	REV.																
C	10363-AP		X01																
						<table border="1" style="width: 100%; border-collapse: collapse;"> <tr> <th style="width: 25%;">SCALE</th> <th style="width: 25%;">NEXT ASSY.</th> <th style="width: 25%;">SHEET</th> <th style="width: 25%;">OF</th> </tr> <tr> <td>1:10</td> <td></td> <td>1</td> <td>3</td> </tr> </table>		SCALE	NEXT ASSY.	SHEET	OF	1:10		1	3				
SCALE	NEXT ASSY.	SHEET	OF																
1:10		1	3																

UTILITY REQUIREMENT NOTES:

1. FLUSH LINES CLEAR BEFORE CONNECTING TO STATION.
2. PLUMBER TO PROVIDE SHUT-OFF VALVES FOR WATER CONNECTIONS.
3. PROVIDE 60-65 PSI WATER PRESSURE AND 4 GALLONS PER MINUTE OF FLOW FOR PROPER OPERATION OF HYDRO ASPIRATOR.
4. ELECTRICIAN TO PROVIDE LIQUID TIGHT JUNCTION BOX AND CONDUIT CONNECTIONS PER LOCAL CODES.


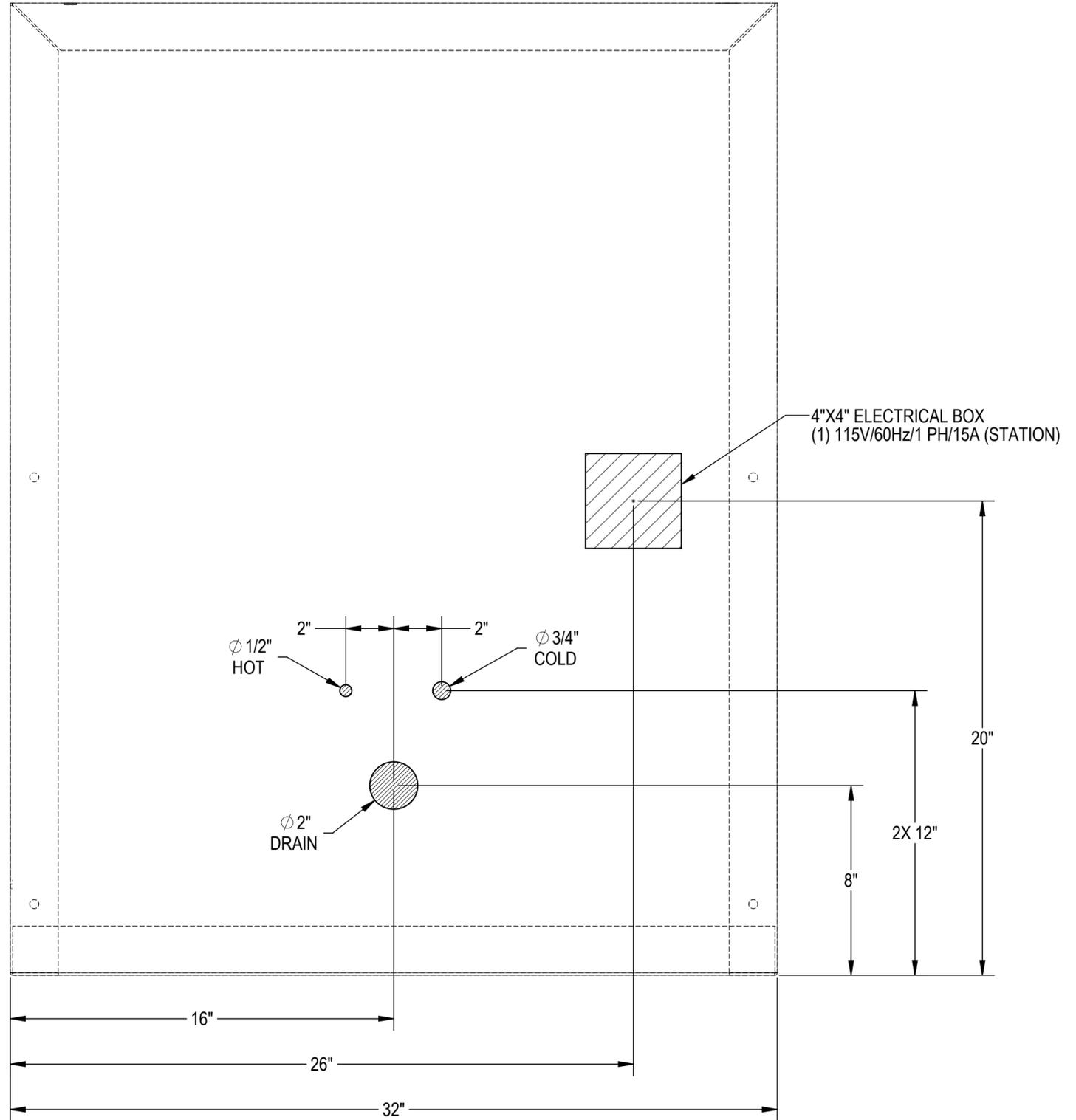


CUSTOMER	APPROVAL		<p style="text-align: center;">PROPRIETARY</p> <p style="text-align: center;">THIS DOCUMENT AND THE DATE DISCLOSED HEREIN AND HEREWITH IS NOT TO BE REPRODUCED, USED, OR OTHERWISE DISCLOSED IN WHOLE OR IN PART TO ANYONE WITHOUT WRITTEN AUTHORIZATION FROM MORTECH MFG.</p>			<p style="font-size: small;">411 North Aerojet Avenue Azusa, CA 91702 tel: (800) 410-0100 (626) 334-1471 fax: (626) 334-1704 www.mortechmfg.com</p>	
	SIGNATURE						
PROJECT	DO NOT SCALE DRAWING		APPROVALS	DATE	SIZE	DWG NO.	REV.
	MATERIAL		CHECKED BY		C		
	304 STAINLESS STEEL		GABRIEL G.	03/10/17			
	FINISH		APPROVED BY		SCALE	NEXT ASSY.	SHEET
	#4 SATIN				1:4		2 OF 3

WALL MOUNTING NOTES:

1. Z-BAR SHIPPED WITHOUT MOUNTING HOLES FOR ADOPTION TO FIELD CONDITIONS. HOLES TO BE DRILLED BY INSTALLER.
2. INSTALL MINIMUM (4) UPPER ANCHORS.
FIRST HOLE: 1" IN FROM SIDE OF Z-BAR.
Z-BAR LENGTH: 27" LONG.
ADDITIONAL ANCHORS CAN BE ADDED AS NEEDED.
3. USE MINIMUM 1/4" TEK SCREWS 2 3/4" LONG. ADJUST LENGTH PER WALL THICKNESS.
4. ANCHORS USED MUST BE FLUSH TO Z-BAR AS MUCH AS POSSIBLE. THEY MUST NOT INTERFERE WITH PLUMBING BEHIND WALL.
5. THE LOWER ANCHORS ARE USED TO KEEP THE BOTTOM OF THE SINK FROM PULLING AWAY FROM THE WALL. THE Z-BAR HOLDS THE WEIGHT OF THE SINK.
6. INSTALLATION IS BASED ON ROUGH-IN MOUNTING DETAIL PROVIDED.

SUPPORT LOCATION FOR EMBALMING STATION ONLY.
ADDITIONAL WALL SUPPORTS MAY BE REQUIRED.

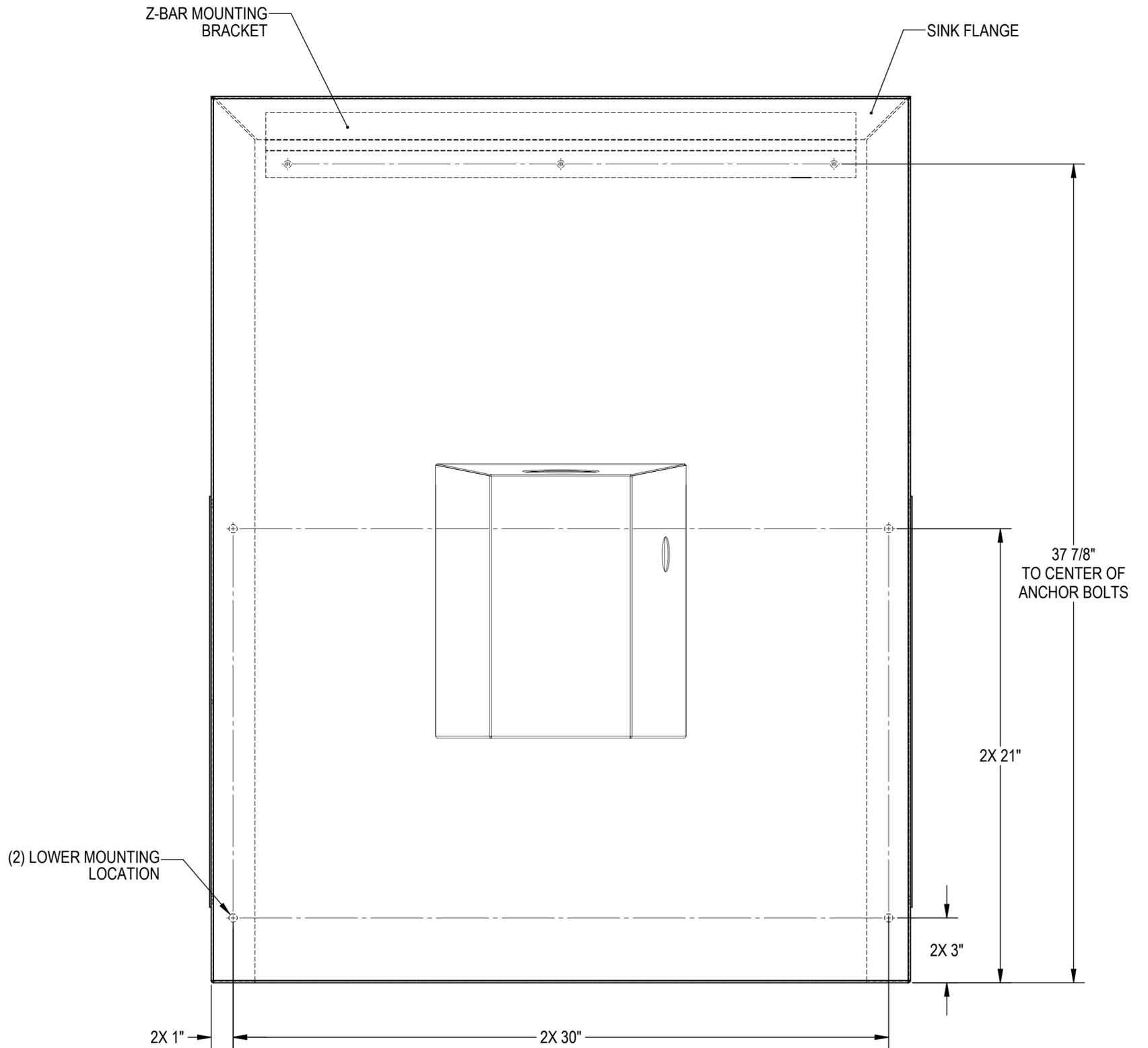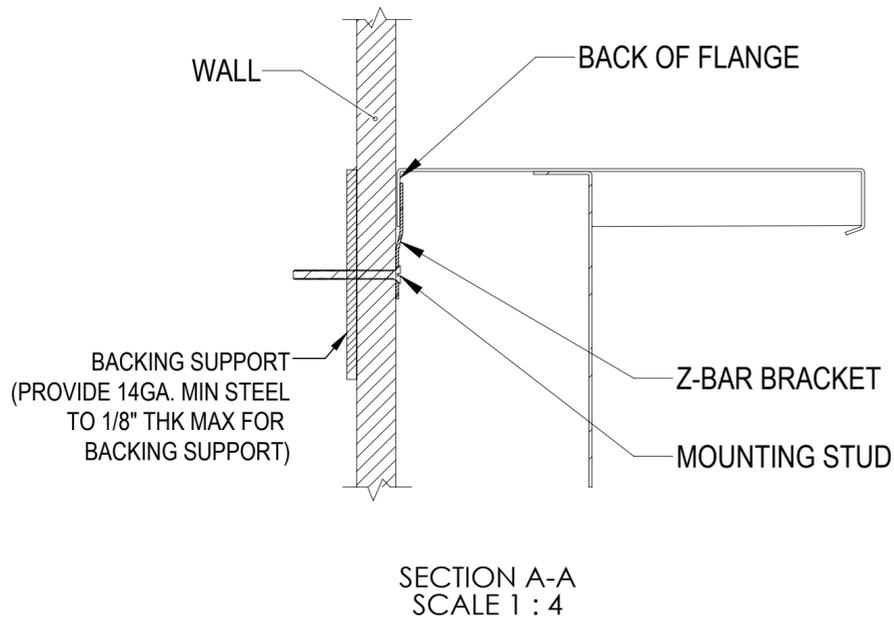

CUSTOMER	APPROVAL	PROPRIETARY			 411 North Aerojet Avenue Azusa, CA 91702 tel: (800) 410-0100 (626) 334-1471 fax: (626) 334-1704 www.mopecmfg.com
		THIS DOCUMENT AND THE DATE DISCLOSED HEREIN AND HEREWITH IS NOT TO BE REPRODUCED, USED, OR OTHERWISE DISCLOSED IN WHOLE OR IN PART TO ANYONE WITHOUT WRITTEN AUTHORIZATION FROM MORTECH MFG.			
PROJECT	SIGNATURE	DATE	DO NOT SCALE DRAWING	APPROVALS	DATE
			MATERIAL	CHECKED BY	REV.
			304 STAINLESS STEEL	GABRIEL G.	03/10/17
			FINISH	APPROVED BY	
			#4 SATIN		
			TITLE MODEL 1036-3 STAINLESS STEEL EMBALMING STATION W/ EMBALMING COLLECTOR		DWG NO. 10363-AP
			SCALE	NEXT ASSY.	REV. X01
			1:4		
			SHEET	OF	
			3	3	