

REV	DESCRIPTION	DATE	ECN #	APPR
A	INITIAL RELEASE	00-XXX-19	XXXXX	MM

LOWERED

RAISED



LA100/OA111 WIRING WORK INSTRUCTIONS SEE: INS-026

Item Number	Quantity	File Name	Title	Material	Revision
1'	6	OA111-001.par	ADAPTER FRAME SPACERS OPTIONAL		A
2'	1	OA111-002.psm	OA111 PANEL MOUNT BRACKET	SHT, SS 11 GA, 304	A
3	1	PE0007.par	PE0007.par		A
4	2	PE0078.par	CABLE GRIP 1/2 NPT		A
5	2	PE0079.par	STRAIN RELIEF LINE CORD .170-.470" Diam, 2" Flex Prot, Black		
6	2	PP0711.par	ORION WASTE PIPE ADAPTER	PROGRESSIVE # 01101518	
7	1	PP0713.par	ORION WASTE PIPE SERRATED FITTING	MCMaster-CARR # 5218K49	
8	1	PP0715.par	ORION WASTE PIPE COUPLER	PROGRESSIVE # 131101501	
9	1	PPI256.par	WASTE DRAIN VALVE, TWIST HANDEL		
10	1	S7952-A.asm	OA110 MOD		A
11	1	S7956-A.asm	OA110 MOD TABLE TOP		A
12	1	S7965-A.psm	J-HOOK	SHT, SS 11 GA, 304	A


<p>THIS DOCUMENT CONTAINS PROPRIETARY INFORMATION OF MOPEC, AND MAY NOT BE COPIED IN WHOLE OR PART, AND NO USE SHALL BE MADE OF IT OTHER THAN FOR WHICH IT HAS BEEN SUBMITTED WITHOUT THE WRITTEN PERMISSION OF A MOPEC MANAGER</p> <p>MOPEC 2020</p>	<p>UNLESS OTHERWISE SPECIFIED ALL DIMENSIONS IN INCHES</p> <p>TOLERANCES</p> <table> <tr> <td>0.0</td> <td>± 0.1 in</td> </tr> <tr> <td>0.00</td> <td>± 0.05 in</td> </tr> <tr> <td>0.000</td> <td>± 0.010 in</td> </tr> <tr> <td>0.0000</td> <td>± 0.0010 in</td> </tr> <tr> <td>ANGLES</td> <td>± 0.5°</td> </tr> </table> <p>BREAK ALL CORNERS 0.063 X 45° ALL FILLET RADII R0.063 (0.125in) INSPECT OVAL DIMENSIONS</p>	0.0	± 0.1 in	0.00	± 0.05 in	0.000	± 0.010 in	0.0000	± 0.0010 in	ANGLES	± 0.5°	<p>MATERIAL</p> <p>MATERIAL USAGE</p> <p>Error: No reference</p> <p>FINISH</p> <p>Error: No reference</p>	<p><b>MOPEC GROUP</b></p> <p>800 Tech Row, Madison Heights, MI 48071</p> <p>DRAWN TO 3RD ANGLE PROJECTION</p>	<p>DRAWN BY: dallmayer DATE: 2/12/2026</p> <p>CHECKED BY: --- DATE: 2/12/2026</p> <p>APPROVED BY: MMcCLAIN DATE: 2/12/2026</p>	<p>PART NUMBER: LA100-A</p> <p>DESCRIPTION: OA111 LARGE NECROPSY TABLE</p> <p>SIZE: A2 FILE NAME: LA100-A.dft</p> <p>SCALE: 1:6.67</p>	<p>DO NOT SCALE IF IN DOUBT ASK</p>	<p>REV: A</p> <p>SHEET 1 OF 4</p>
		0.0	± 0.1 in														
0.00	± 0.05 in																
0.000	± 0.010 in																
0.0000	± 0.0010 in																
ANGLES	± 0.5°																
<p>800 Tech Row, Madison Heights, MI 48071</p>																	

\*LA100-A.dft\*


REV	DESCRIPTION	DATE	ECN #	APPR
A	INITIAL RELEASE	00-XXX-19	XXXXX	MM

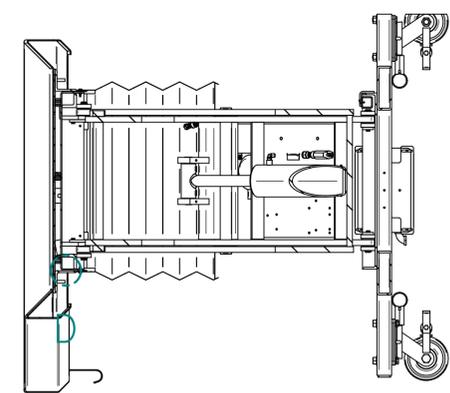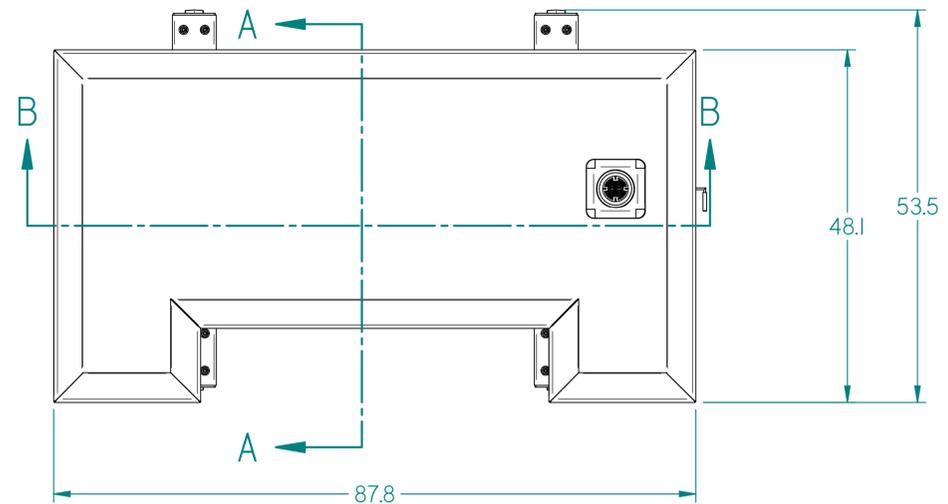


SECTION A-A

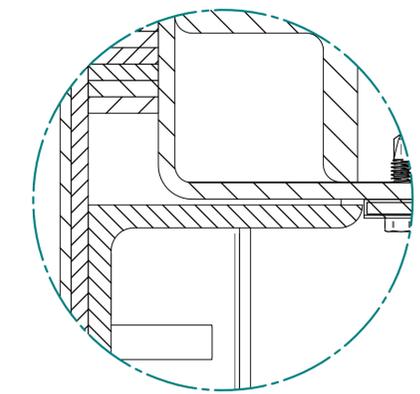



DETAIL D

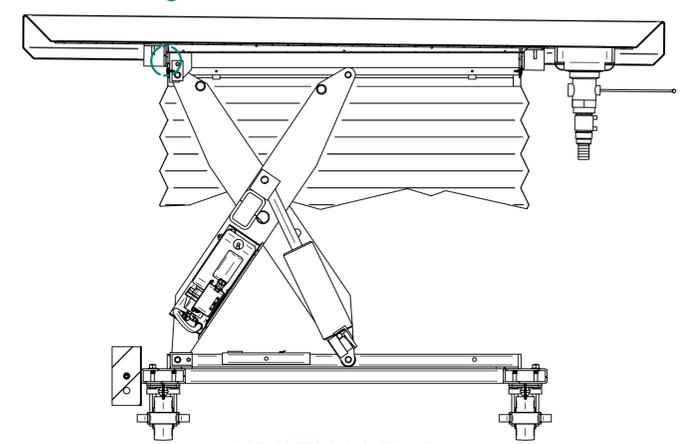

SECTION B-B

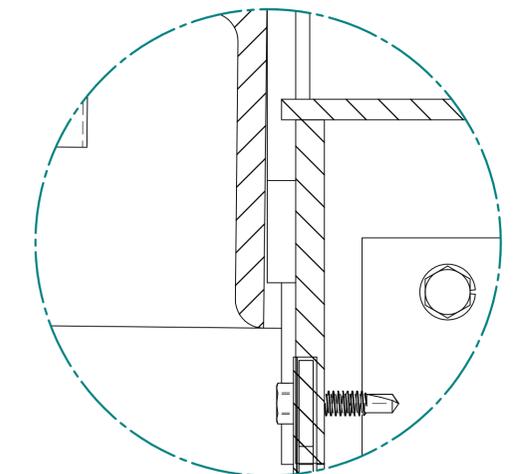



DETAIL C

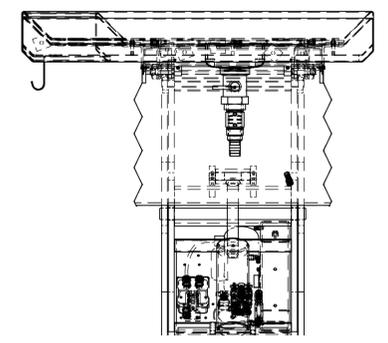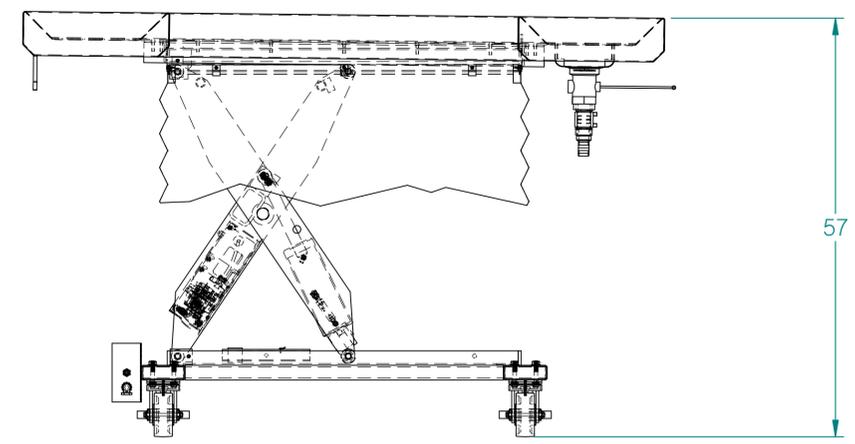


<p>THIS DOCUMENT CONTAINS PROPRIETARY INFORMATION OF MOPEC, AND MAY NOT BE COPIED IN WHOLE OR PART, AND NO USE SHALL BE MADE OF IT OTHER THAN FOR WHICH IT HAS BEEN SUBMITTED WITHOUT THE WRITTEN PERMISSION OF A MOPEC MANAGER</p> <p>MOPEC 2020</p>	<p>UNLESS OTHERWISE SPECIFIED ALL DIMENSIONS IN INCHES</p> <p>TOLERANCES</p> <table> <tr><td>0.0</td><td>± 0.1 in</td></tr> <tr><td>0.00</td><td>± 0.05 in</td></tr> <tr><td>0.000</td><td>± 0.010 in</td></tr> <tr><td>0.0000</td><td>± 0.0010 in</td></tr> <tr><td>ANGLES</td><td>± 0.5°</td></tr> </table> <p>BREAK ALL CORNERS 0.063 X 45° ALL FILLET RADII R0.063 (0.125in) INSPECT OVAL DIMENSIONS</p>	0.0	± 0.1 in	0.00	± 0.05 in	0.000	± 0.010 in	0.0000	± 0.0010 in	ANGLES	± 0.5°	<p>MATERIAL</p> <p>MATERIAL USAGE</p> <p>Error: No reference</p> <p>FINISH</p> <p>Error: No reference</p>	<p><b>MOPEC GROUP</b></p> <p>800 Tech Row, Madison Heights, MI 48071</p> <p>DRAWN TO 3RD ANGLE PROJECTION</p>	<p>DRAWN BY: dallmayer</p> <p>DATE: 2/12/2026</p> <p>CHECKED BY: ---</p> <p>DATE: 2/12/2026</p> <p>APPROVED BY: M.McCLAIN</p> <p>DATE: 2/12/2026</p>	<p>PART NUMBER: LA100-A</p> <p>DESCRIPTION: OAH LARGE NECROPSY TABLE</p> <p>SIZE: A2</p> <p>SCALE: 1:8</p> <p>DO NOT SCALE IF IN DOUBT ASK</p>	<p>FILE NAME: LA100-A.dft</p> <p>REV: A</p> <p>SHEET 2 OF 4</p>
	0.0	± 0.1 in														
0.00	± 0.05 in															
0.000	± 0.010 in															
0.0000	± 0.0010 in															
ANGLES	± 0.5°															
<p>UNLESS OTHERWISE SPECIFIED ALL DIMENSIONS IN INCHES</p>		<p>800 Tech Row, Madison Heights, MI 48071</p>		<p>LA100-A</p>												

\*LA100-A.dft\*


REV	DESCRIPTION	DATE	ECN #	APPR
A	INITIAL RELEASE	00-XXX-19	XXXXX	MM

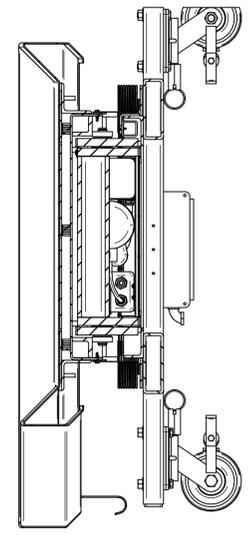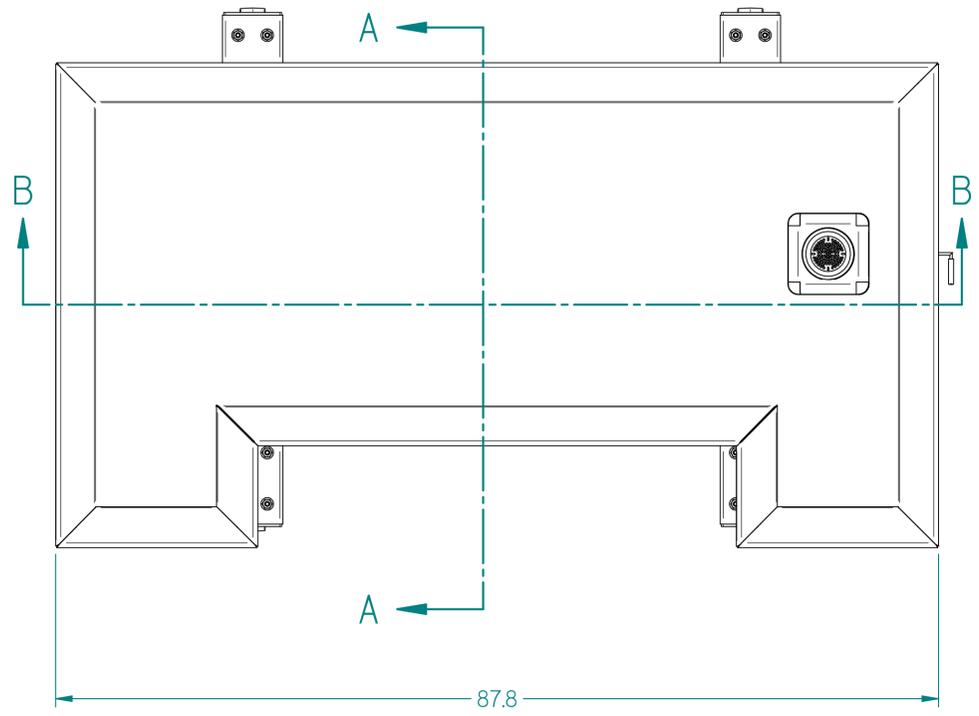

SECTION A-A

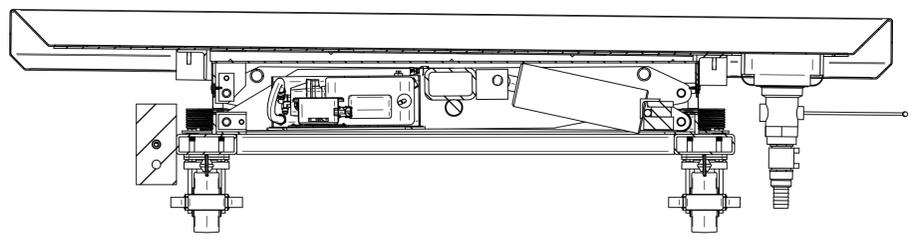


SECTION B-B

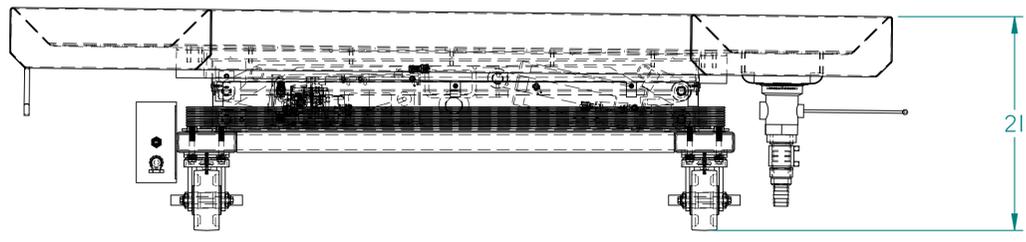

<p>THIS DOCUMENT CONTAINS PROPRIETARY INFORMATION OF MOPEC, AND MAY NOT BE COPIED IN WHOLE OR PART, AND NO USE SHALL BE MADE OF IT OTHER THAN FOR WHICH IT HAS BEEN SUBMITTED WITHOUT THE WRITTEN PERMISSION OF A MOPEC MANAGER</p> <p>MOPEC 2020</p>	<p>UNLESS OTHERWISE SPECIFIED ALL DIMENSIONS IN INCHES</p> <p>TOLERANCES</p> <table> <tr><td>0.0</td><td>± 0.1 in</td></tr> <tr><td>0.00</td><td>± 0.05 in</td></tr> <tr><td>0.000</td><td>± 0.010 in</td></tr> <tr><td>0.0000</td><td>± 0.0010 in</td></tr> </table> <p>ANGLES ±0.5°</p> <p>BREAK ALL CORNERS 0.063 X 45°</p> <p>ALL FILLET RADII R0.063 (0.125in)</p> <p>INSPECT OVAL DIMENSIONS</p>	0.0	± 0.1 in	0.00	± 0.05 in	0.000	± 0.010 in	0.0000	± 0.0010 in	<p>MATERIAL</p> <p>MATERIAL USAGE</p> <p>Error: No reference</p> <p>FINISH</p> <p>Error: No reference</p>	<p><b>MOPEC GROUP</b></p> <p>800 Tech Row, Madison Heights, MI 48071</p> <p>DRAWN BY: dallmayer DATE: 2/12/2026</p> <p>CHECKED BY: --- DATE: ---/---/2026</p> <p>APPROVED BY: M.McCLAIN DATE: 2/12/2026</p>	<p>PART NUMBER: LA100-A</p> <p>DESCRIPTION: OAH LARGE NECROPSY TABLE</p> <p>SIZE: A2 FILE NAME: LA100-A.dft</p> <p>SCALE: 1:8</p> <p>DO NOT SCALE IF IN DOUBT ASK</p>	<p>REV: A</p> <p>SHEET 3 OF 4</p>
	0.0	± 0.1 in											
0.00	± 0.05 in												
0.000	± 0.010 in												
0.0000	± 0.0010 in												
<p>DRAWN TO 3RD ANGLE PROJECTION</p>		<p>800 Tech Row, Madison Heights, MI 48071</p>		<p>800 Tech Row, Madison Heights, MI 48071</p>									


\*LA100-A.dft\*

REV	DESCRIPTION	DATE	ECN #	APPR
A	INITIAL RELEASE	00-XXX-19	XXXXX	MM

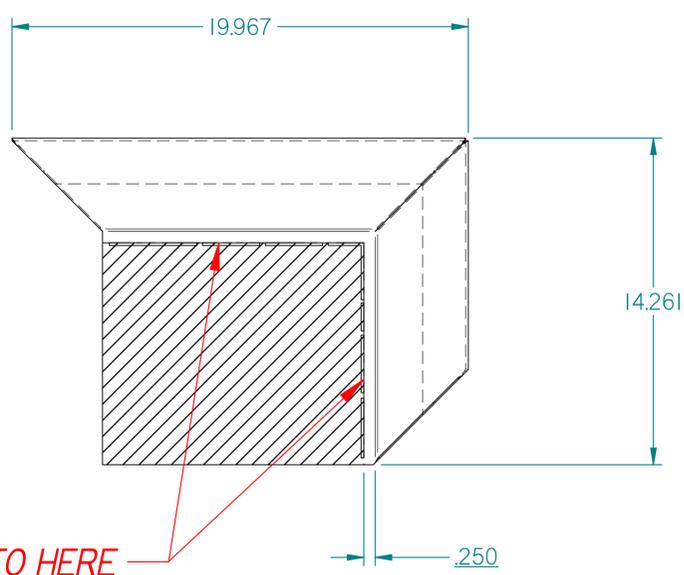

CUT BACK MATERIAL TO HERE AFTER FORMING

CUT BACK MATERIAL AFTER FORMING TO CLEAR I36253-I-121-A BACKING PLATE

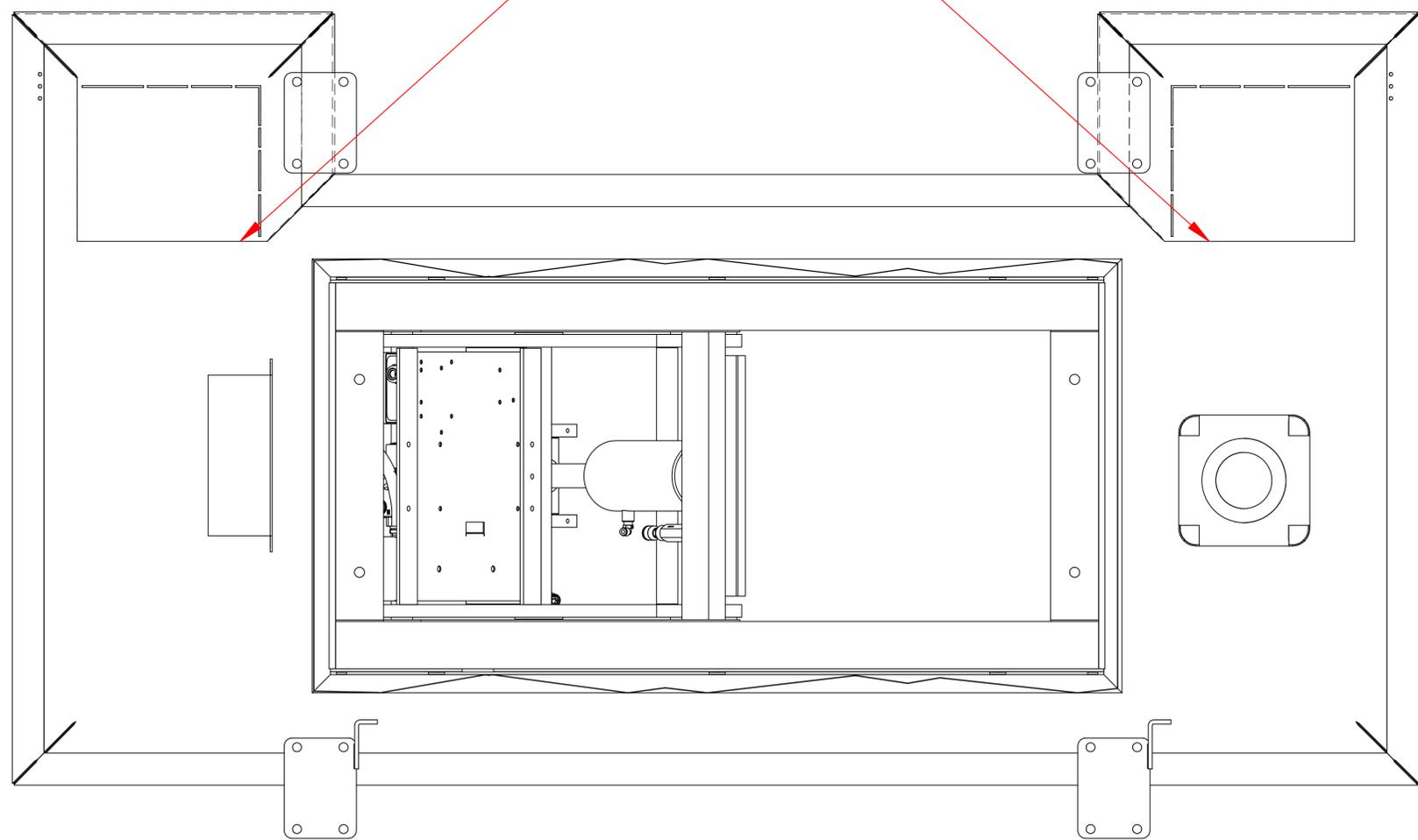

<p>THIS DOCUMENT CONTAINS PROPRIETARY INFORMATION OF MOPEC, AND MAY NOT BE COPIED IN WHOLE OR PART, AND NO USE SHALL BE MADE OF IT OTHER THAN FOR WHICH IT HAS BEEN SUBMITTED WITHOUT THE WRITTEN PERMISSION OF A MOPEC MANAGER</p> <p>MOPEC 2020</p>	<p>UNLESS OTHERWISE SPECIFIED ALL DIMENSIONS IN INCHES</p> <p>TOLERANCES</p> <table> <tr><td>0.0</td><td>± 0.1 in</td></tr> <tr><td>0.00</td><td>± 0.05 in</td></tr> <tr><td>0.000</td><td>± 0.010 in</td></tr> <tr><td>0.0000</td><td>± 0.0010 in</td></tr> <tr><td>ANGLES</td><td>± 0.5°</td></tr> </table> <p>BREAK ALL CORNERS 0.063 X 45° ALL FILLET RADII R0.063 (0.125in) INSPECT OVAL DIMENSIONS</p>	0.0	± 0.1 in	0.00	± 0.05 in	0.000	± 0.010 in	0.0000	± 0.0010 in	ANGLES	± 0.5°	<p>MATERIAL</p> <p>MATERIAL USAGE</p> <p>Error: No reference</p> <p>FINISH</p> <p>Error: No reference</p>	<p><b>MOPEC GROUP</b></p> <p>800 Tech Row, Madison Heights, MI 48071</p> <p>DRAWN TO 3RD ANGLE PROJECTION</p>	<p>DRAWN BY: dallmayer DATE: 2/12/2026</p> <p>CHECKED BY: --- DATE: ---/---/2026</p> <p>APPROVED BY: MMcCLAIN DATE: 2/12/2026</p>	<p>PART NUMBER: LA100-A</p> <p>DESCRIPTION: OAH LARGE NECROPSY TABLE</p> <p>SIZE: A2</p> <p>SCALE: 1:1</p> <p>DO NOT SCALE IF IN DOUBT ASK</p>	<p>REV: A</p> <p>SHEET 4 OF 4</p>
	0.0	± 0.1 in														
0.00	± 0.05 in															
0.000	± 0.010 in															
0.0000	± 0.0010 in															
ANGLES	± 0.5°															
<p>FILE NAME: LA100-A.dft</p>		<p>800 Tech Row, Madison Heights, MI 48071</p>		<p>LA100-A</p>												

\*LA100-A.dft\*