


REV	DESCRIPTION	DATE	ECN #	APPR
B	PARTIAL, STAINLESS PRODUCT RELEASE	10-MAR-15	15814	KLB

SECTION A-A

Item #	Qty.	File Name	Title	Rev.
1'	2	PE0015-A.par	COVER RECEPTACLE GFCI, 1-Gang, Stay-Open Lid	A
2'	1	PE0016.par	COVER PLUG BLANK, 1-GANG, SS W/SCREWS & GASKET	
3'	3	PE0309-A.par	BOX ELEC. OUTLET-WATERPROOF 2-1/2" DEEP	A
4'	1	PE0438.par	WATER PROOF 20AMP C/B SWITCH (NON-DISPOSAL)	A
5'	2	PE0510.par	RECEPTACLE GFCI, 20 AMP, 125V, Str Blade, Gray, Hospital	A
6'	2	PP0034.par	STAINLESS STEEL DRAIN STRAINER - 1-1/2" NPT	
7'	2	PP0058.par	FITTING - BRASS TAIL DRAIN TUBE - 1-1/2" x 4"L	A
8	2	PP0065.par	FITTING - SPOUT END - B SPOUT	
9'	1	PP0140.asm	HOSEASSY1.asm	
10'	2	PP0199.par	FITTING BRASS B/F PREVENTER 3/8" FNPT	A
11	1	PP0238.par	FISHER 3700 SINGLE CONTROL VALVE	
12	2	PP0295.par	FIXTURE - CHROME SPOUT - 6" INLET	A
13	4	PP0400.par	FIXTURE HANDLES	
14	1	PP0507.par	FIXTURE - 4" CENTER CHICAGO	A
15	1	PP0520.par	FIXTURE - FAUCET W/O SPOUT	A
16'	1	SI230-C.psm	J-HOOK	C
17	3	SI839-B.psm	COMMON OA TUB BAFFLE	B
18	3	SI840-B.psm	COMMON OA GRID PLATE	A
19	1	SI841-B.asm	OA PLENUM WLDT	B
20	1	SI849-B.asm	OA SERIES DAMPENR ASM	B
21	1	S6072-B.psm	UTILITY COVER	B
22	1	S6073-A.asm	OA800 MAJOR WLDT	A

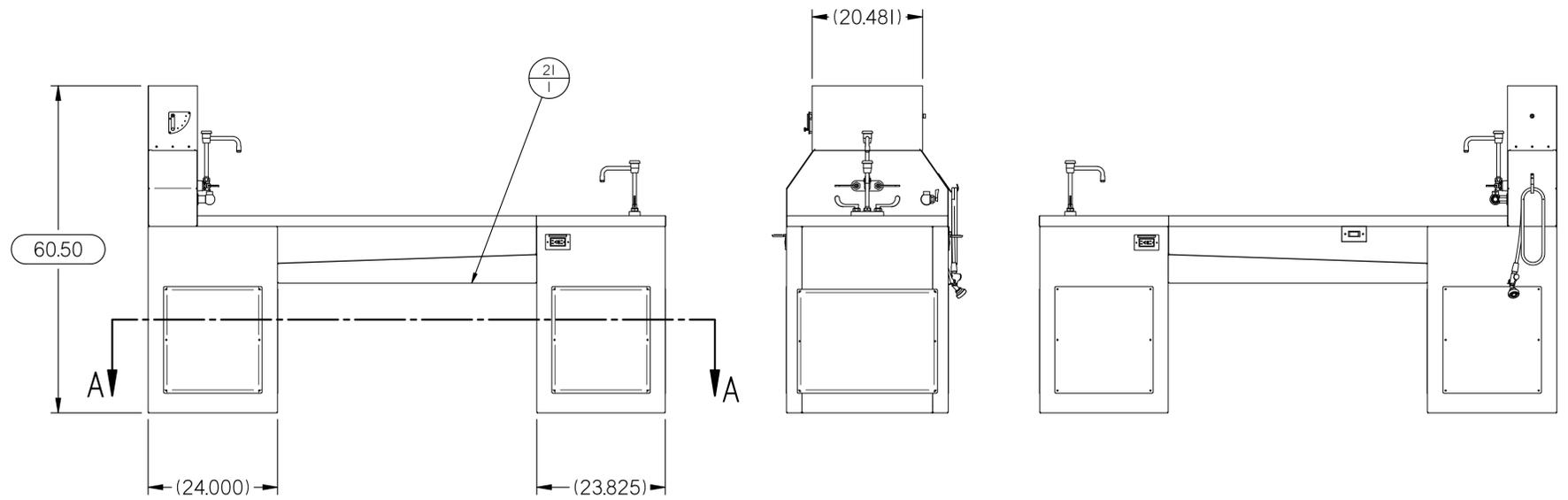

MOPEC LEGAL NOTE THIS DOCUMENT CONTAINS PROPRIETARY INFORMATION OF MOPEC, AND MAY NOT BE COPIED IN WHOLE OR PART, AND NO USE SHALL BE MADE OF IT OTHER THAN FOR WHICH IT HAS BEEN SUBMITTED WITHOUT THE WRITTEN PERMISSION OF A MOPEC MANAGER MOPEC 2025	UNLESS OTHERWISE SPECIFIED ALL DIMENSIONS IN INCHES TOLERANCES 0.0 ± 0.1 in 0.00 ± 0.05 in 0.000 ± 0.010 in 0.0000 ± 0.0010 in ANGLES ±0.5° BREAK ALL CORNERS 0.063 X 45° ALL FILLET RADII R0.063 /0.125in INSPECT OVAL DIMENSIONS	 	DRAWN BY: kbuttery DATE: 1/30/2020	PART NUMBER: OA800-B DESCRIPTION: OA800 NECROPSY TABLE
			CHECKED BY: DATE: 1/30/2020	FILE NAME: OA800-B.dft SCALE: 1:15 DO NOT SCALE IF IN DOUBT ASK
DRAWN TO 3RD ANGLE PROJECTION			APPROVED BY: M.McCLAIN DATE: 1/30/2020	SHEET 1 OF 2

OA800-B.dft

