


REV	DESCRIPTION	DATE	ECN #	APPR
A	INITIAL RELEASE	00-XXX-19	XXXXX	MM
B	PARTIAL, STAINLESS PRODUCT RELEASE	10-MAR-15	15814	KLB
C	STANDARDIZED	27-APR-21	1045	MM
D	LINED UP HOLES	13-MAY-21	1047	MM




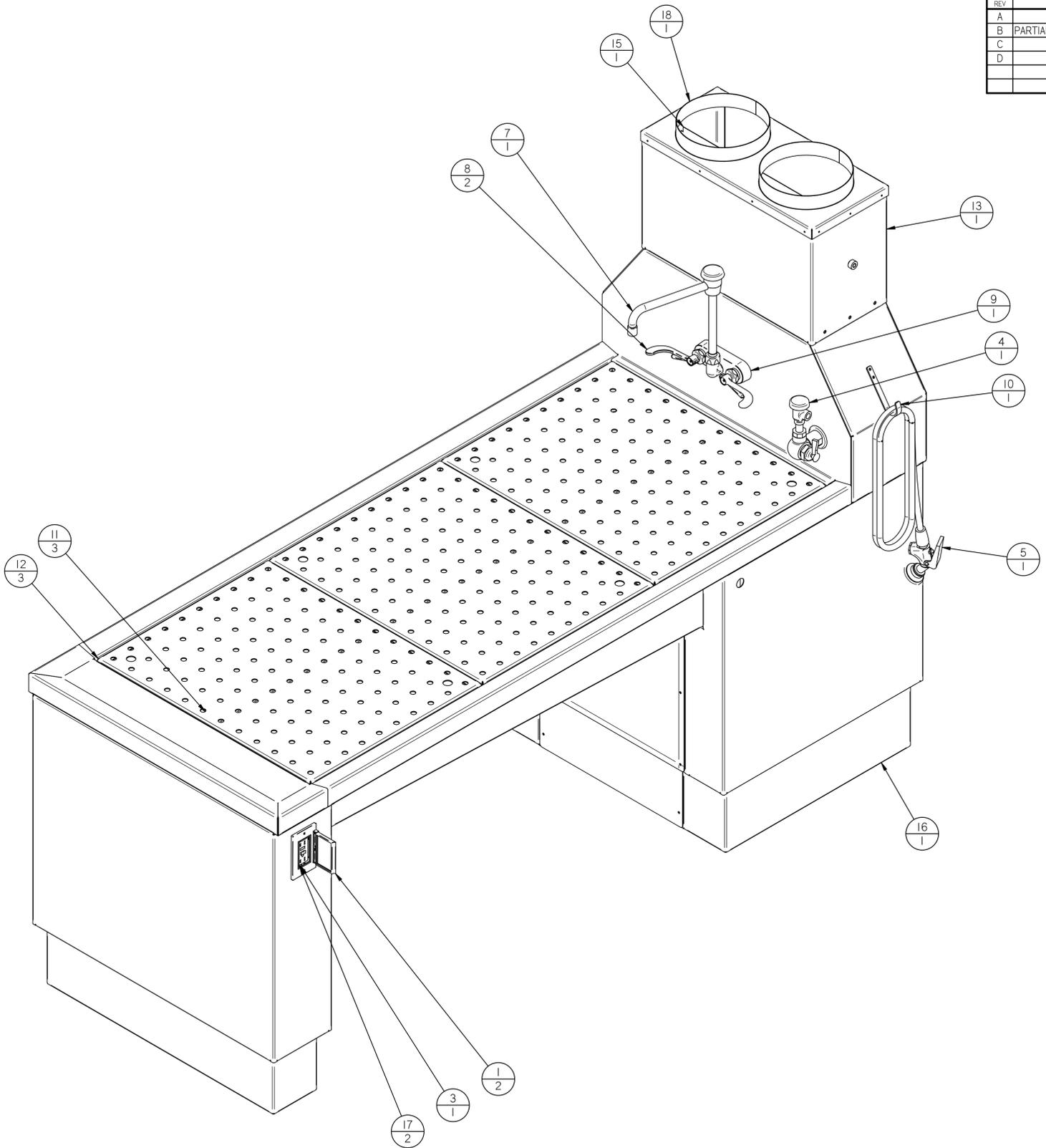


Item Number	Quantity	File Name	Title	Revision
1	2	PE0015-A.par	COVER RECEPTACLE GFCI, 1-Gang, Stay-Open Lid	A
2	1	PE0249.par	RECEPTACLE, 20 Amp, 125V Duplex, Hospital Grade, Gray	A
3	1	PE0510.par	RECEPTACLE GFCI, 20 AMP, 125V, Str Blade, Gray, Hospital	A
4	1	PP0095.par	FIXTURE - BRASS VAC. BREAKER 90 DEG - 1/4" FNPT OUTLET	A
5	1	PP0140.asm	HOSEASSY1.asm	
6	1	PP0238.par	FISHER 3700 SINGLE CONTROL VALVE	
7	1	PP0294.par	FIXTURE - CHROME SPOUT - 8" INLET	A
8	2	PP0400.par	FIXTURE HANDLES	
9	1	PP0520.asm	PP0520.asm	A
10	1	SI230-C.psm	J-HOOK	C
11	3	SI839-B.psm	COMMON OA TUB BAFFLE	B
12	3	SI840-B.psm	COMMON OA GRID PLATE	A
13	1	SI841-B.asm	OA PLENUM WLDT	B
14	1	SI845-C.psm	OA UTILITY COVER - SHORT	C
15	1	SI849-B.asm	OA SERIES DAMPENER ASM	B
16	1	S2200-C.asm	OA900 MAJOR WLDT	C
17	2	S2960-A.psm	DUPLEX BOX ELIMINATOR PLATE	A
18	1	S6077-A.asm	OA900 DUCT MANIFOLD WLDT	A



THIS DOCUMENT CONTAINS PROPRIETARY INFORMATION OF MOPEC, AND MAY NOT BE COPIED IN WHOLE OR PART, AND NO USE SHALL BE MADE OF IT OTHER THAN FOR WHICH IT HAS BEEN SUBMITTED WITHOUT THE WRITTEN PERMISSION OF A MOPEC MANAGER MOPEC 2025	UNLESS OTHERWISE SPECIFIED ALL DIMENSIONS IN INCHES TOLERANCES 0.0 ± 0.1 in 0.00 ± 0.05 in 0.000 ± 0.010 in 0.0000 ± 0.0010 in ANGLES ±0.5° BREAK ALL CORNERS 0.063 X 45° ALL FILLET RADII R0.063 / 0.125 in INSPECT OVAL DIMENSIONS	 MOPEC GROUP 800 Tech Row, Madison Heights, MI 48071	DRAWN BY: jmazza DATE: 5/13/2021	PART NUMBER: OA900-D
			CHECKED BY: --- DATE: 5/13/2021	DESCRIPTION: TRIMMING TABLE - TOTAL DRAFT ELEVATING
DRAWN TO 3RD ANGLE PROJECTION			APPROVED BY: M.McCLAIN DATE: 5/13/2021	SIZE: A2 FILE NAME: OA900-D.dft SCALE: 1:10 DO NOT SCALE IF IN DOUBT ASK

\*OA900-D.dft\*

REV	DESCRIPTION	DATE	ECN #	APPR
A	INITIAL RELEASE	00-XXX-19	XXXXX	MM
B	PARTIAL, STAINLESS PRODUCT RELEASE	10-MAR-15	15814	KLB
C	STANDARDIZED	27-APR-21	1045	MM
D	LINED UP HOLES	13-MAY-21	1047	MM

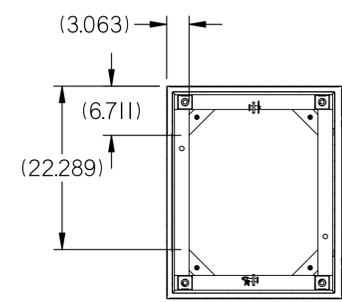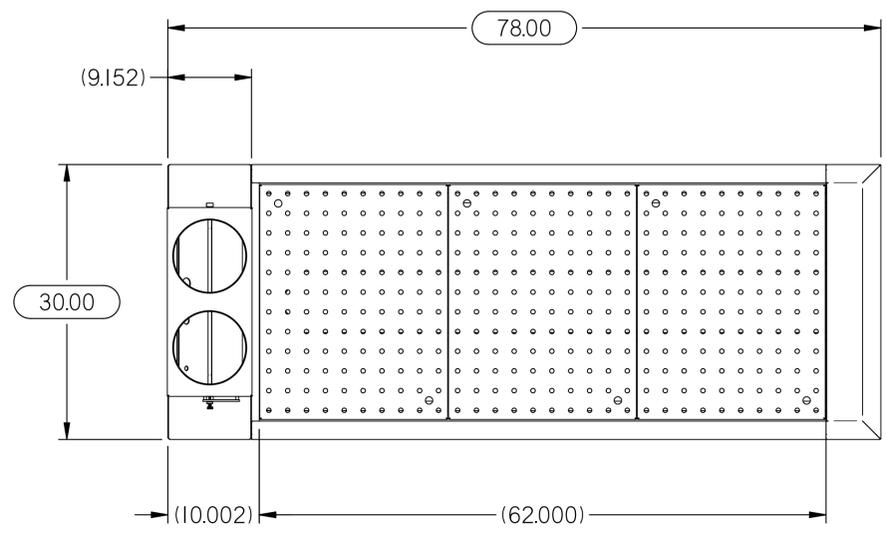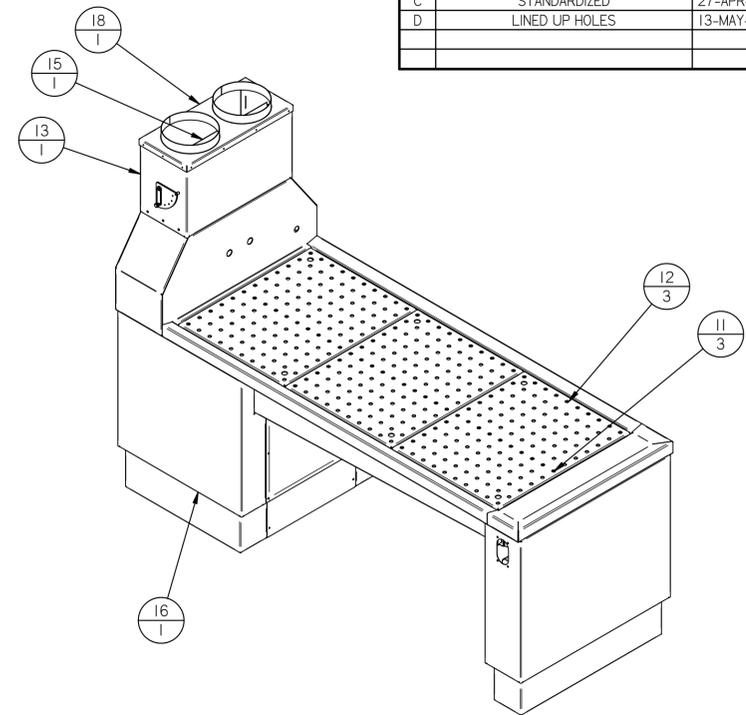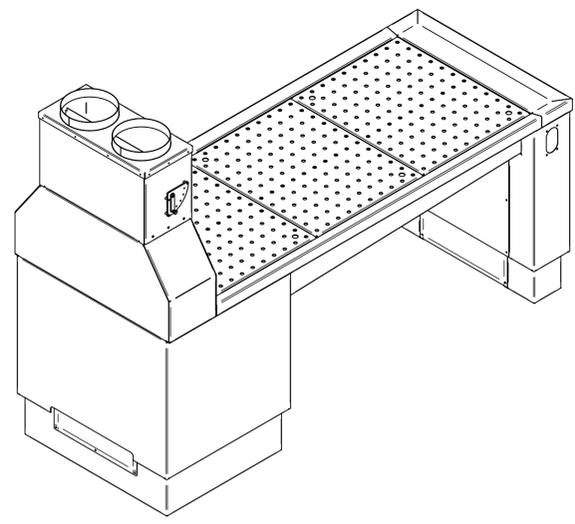

SECTION A-A

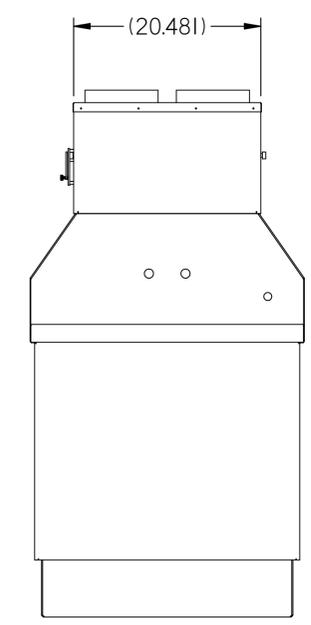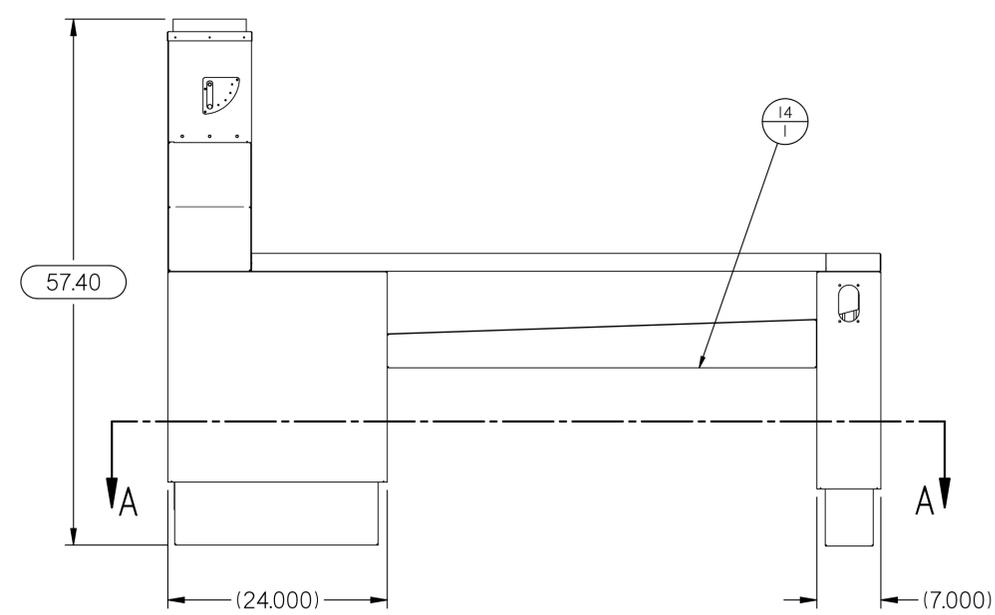

Item Number	Quantity	File Name	Title	Revision
1	2	PE0015-A.par	COVER RECEPTACLE GFCI, I-Gang, Stay-Open Lid	A
2	1	PE0249.par	RECEPTACLE, 20 Amp, 125V Duplex, Hospital Grade, Gray	A
3	1	PE0510.par	RECEPTACLE GFCI, 20 AMP, 125V, Str Blade, Gray, Hospital	A
4	1	PP0095.par	FIXTURE - BRASS VAC. BREAKER 90 DEG - 1/4" FNPT OUTLET	A
5	1	PP0140.asm	HOSEASSY1.asm	
6	1	PP0238.par	FISHER 3700 SINGLE CONTROL VALVE	
7	1	PP0294.par	FIXTURE - CHROME SPOUT - 8" INLET	A
8	2	PP0400.par	FIXTURE HANDLES	
9	1	PP0520.asm	PP0520.asm	A
10	1	SI230-C.psm	J-HOOK	C
11	3	SI839-B.psm	COMMON OA TUB BAFFLE	B
12	3	SI840-B.psm	COMMON OA GRID PLATE	A
13	1	SI841-B.asm	OA PLENUM WLDT	B
14	1	SI845-C.psm	OA UTILITY COVER - SHORT	C
15	1	SI849-B.asm	OA SERIES DAMPENR ASM	B
16	1	S2200-C.asm	OA900 MAJOR WLDT	C
17	2	S2960-A.psm	DUPLEX BOX ELIMINATOR PLATE	A
18	1	S6077-A.asm	OA900 DUCT MANIFOLD WLDT	A

THIS DOCUMENT CONTAINS PROPRIETARY INFORMATION OF MOPEC, AND MAY NOT BE COPIED IN WHOLE OR PART, AND NO USE SHALL BE MADE OF IT OTHER THAN FOR WHICH IT HAS BEEN SUBMITTED WITHOUT THE WRITTEN PERMISSION OF A MOPEC MANAGER MOPEC 2025	UNLESS OTHERWISE SPECIFIED ALL DIMENSIONS IN INCHES TOLERANCES 0.0 ± 0.1 in 0.00 ± 0.05 in 0.000 ± 0.010 in 0.0000 ± 0.0010 in ANGLES ±0.5° BREAK ALL CORNERS 0.063 X 45° ALL FILLET RADII R0.063 / 0.125 in INSPECT OVAL DIMENSIONS	<b>MOPEC GROUP</b> 800 Tech Row, Madison Heights, MI 48071	DRAWN BY: jmazza DATE: 5/13/2021 CHECKED BY: --- DATE: 5/13/2021 APPROVED BY: M.McCLAIN DATE: 5/13/2021	PART NUMBER: OA900-D DESCRIPTION: TRIMMING TABLE - TOTAL DRAFT ELEVATING SIZE: A2 FILE NAME: OA900-D.dft SCALE: 1:10 DO NOT SCALE IF IN DOUBT ASK	SHEET 2 OF 2
		DRAWN TO 3RD ANGLE PROJECTION	REVISION: D		

\*OA900-D.dft\*